
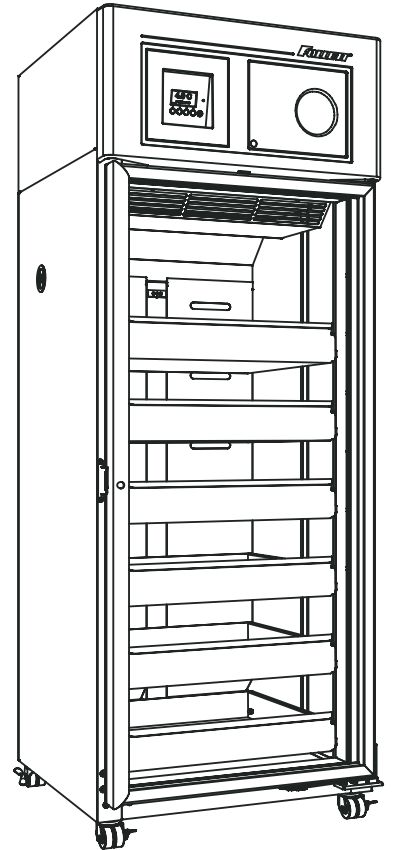
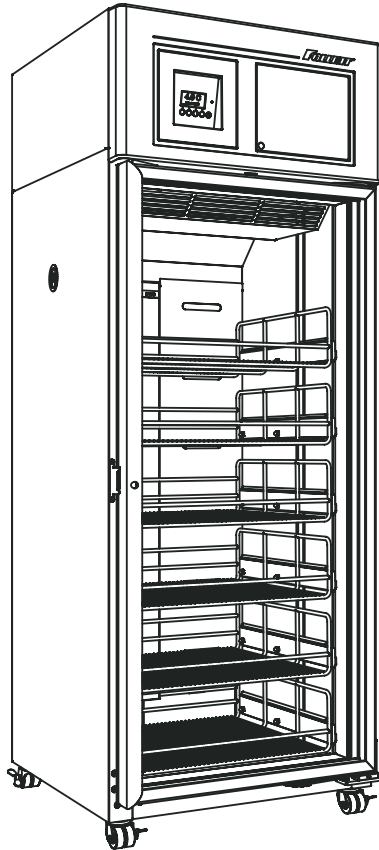


REF 20/25 – LB
REF 20/25 – PH
REF 20/25 – BB

Order parts online
www.follettice.com

Installation, Operation and Service Manual

Serial numbers J62763 and below

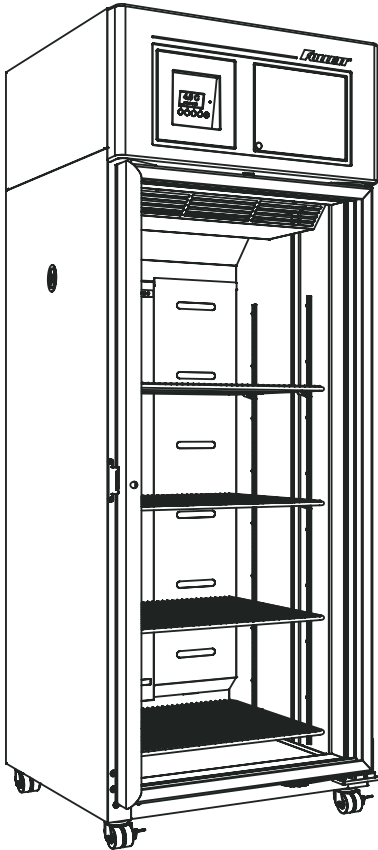

**Following installation, please forward this manual
to the appropriate operations person.**

Contents

Welcome to Follett	3	Cleaning	20
Before you begin	3	Interior cleaning	20
Ambients	4	Exterior cleaning	20
Clearances	4	Annual cleaning	21
Electrical specifications	4	Service	22
Refrigeration specifications	4	Refrigeration system	22
Detail drawing	4	Defrosting	22
Installation	5	Controller operation	23
Adjust self-closing door (if necessary)	5	Temperature sensor readings	23
Install shelves (if equipped) and power up	5	Heated door (option) – changing cycle	24
Fill product bottle	6	Controller hot key	27
Remote alarm connections (if desired)	7	Hot key upload	27
Dry contacts	7	Hot key download	28
RS-485 port	7	Troubleshooting	34
Operation	8	Before calling for service	34
Quick setup checklist	8	Appendix	35
Interior LED lights	8	Control key functions	35
Set time & date	8	Menu Navigation table	35
Change temperature display to show degrees Fahrenheit (if desired)	9	Replacement parts	36
High and low alarm setup	10	Unit cooler	36
Alarms	12	Cabinet	37
Alarm muting	12	Facade	38
Alarm volume selection	13	Door assembly	39
Alarm silence	13		
Password PIN security (default is 3843)	14		
To turn password on/off	14		
To make a new password PIN	15		
Min/max temperature logging	16		
Probe calibration	17		
Chart recorder set up (if so equipped)	18		
Chart recorder calibration	18		
Changing the chart paper	18		
Replacing the pen	18		
Removing drawers	19		
Removing slides	19		


Welcome to Follett

Follett equipment enjoys a well-deserved reputation for excellent performance, long-term reliability and outstanding after-the-sale support. To ensure that this product delivers that same degree of service, we ask that you take a moment to review this manual before beginning the installation. Should you have any questions or require technical help at any point, please call our technical service group at (877) 612-5086 or +1 (610) 252-7301.

Before you begin

After uncrating and removing all packing material, inspect the equipment for concealed shipping damage. If damage is found, notify the shipper immediately and contact Follett Corporation so that we can help in the filing of a claim, if necessary.

Check your paperwork to determine which configuration you have. Follett configuration numbers are designed to provide information about the type of refrigerator you are receiving. Following is an explanation of the different item numbers.

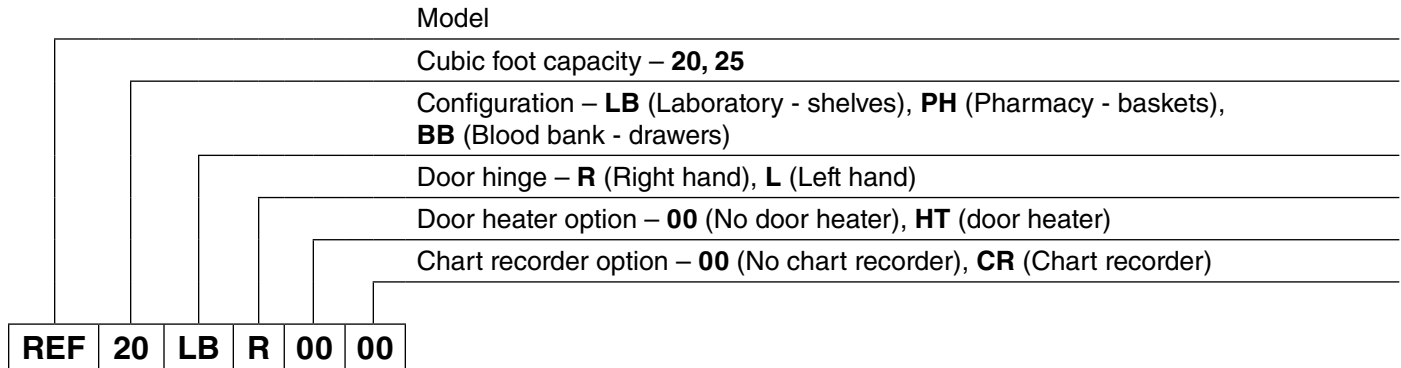

Specifications						
	REF20-LB	REF25-LB	REF20-PH	REF25-PH	REF20-BB	REF25-BB
Capacity	19.7 cu ft (558 L)	24.6 cu ft (697 L)	19.7 cu ft (558 L)	24.6 cu ft (697 L)	19.7 cu ft (558 L)/ 280 bags (450ml)	24.6 cu ft (697 L)/ 360 bags (450ml)
Storage system	(4) epoxy-coated shelves	(4) epoxy-coated shelves	(6) epoxy-coated baskets	(6) epoxy-coated baskets	(6) stainless steel drawers	(6) stainless steel drawers
Exterior width	29.75" (76 cm)	29.75" (76 cm)	29.75" (76 cm)	29.75" (76 cm)	29.75" (76 cm)	29.75" (76 cm)
Exterior depth	29" (74 cm)	35" (89 cm)	29" (74 cm)	35" (89 cm)	29" (74 cm)	35" (89 cm)
Exterior depth with handles	30.5" (78 cm)	36.5" (93 cm)	30.5" (78 cm)	36.5" (93 cm)	30.5" (78 cm)	36.5" (93 cm)
Exterior height with casters	79.5" (202 cm)	79.5" (202 cm)	79.5" (202 cm)	79.5" (202 cm)	79.5" (202 cm)	79.5" (202 cm)
Interior dimensions (w x h x d)	24" x 22.5" x 56" (61 cm x 58 cm x 143 cm)	24" x 28.5" x 56" (61 cm x 73 cm x 143 cm)	24" x 22.5" x 56" (61 cm x 58 cm x 143 cm)	24" x 28.5" x 56" (61 cm x 73 cm x 143 cm)	24" x 22.5" x 56" (61 cm x 58 cm x 143 cm)	24" x 28.5" x 56" (61 cm x 73 cm x 143 cm)
Crated weight	395 (180 kg)	430 (196 kg)	450 (205 kg)	485 (220 kg)	475 (216 kg)	525 (239 kg)
Max. heat rejection	950 BTU/hr	1050 BTU/hr	950 BTU/hr	1050 BTU/hr	950 BTU/hr	1050 BTU/hr

Standard features – all models	
Door	Dual pane, low-E glass, condensation free to 80 F (27 C) air/60% RH
Door handle	ADA-compliant, full length handle with integral lock
Interior/exterior material	Heavy-duty, corrosion resistant stainless steel
Casters	(4) dual-wheel swiveling casters with toe locks
Insulation	2.75" (7 cm) thick, CFC-free foam insulation throughout
Lighting	(2) full length side LED lights
Air circulation system	Ducted air through 12 back plenum openings with front face return
Temperature probe	Stainless steel RTD (resistance temperature detectors) top, bottom probe standard on - BB models, available as option on -LB, -PH models, immersible with panel quick-disconnect

Ambients

Intended for indoor use where ambient temperature, measured at the condenser inlet, does not exceed 80 F (27 C) and ambient relative humidity does not exceed 60% RH. A heated door is recommended for higher humidity ambients.

Clearances

The top of the refrigerator must have 10 inches (25.4 cm) of clear space above the refrigerator to ensure proper ventilation of the refrigeration system.

The back of the refrigerator should have 1 inch (2.5 cm) of clearance to allow for power cord clearance on unit and an additional 1 inch (2.5 cm) if the unit is to be plugged in directly behind the refrigerator.

Electrical specifications

115V, 60Hz, 1 phase

Running load amps: 8.4 amps without heated door; max 9.7 amps with door heater energized 100% (adjustable).

Full load amps (90% maximum continuous current per UL 471 at 40 C): 10.5 amps


Minimum circuit ampacity: 15 amp dedicated circuit

Maximum size of branch circuit overcurrent device: 15 amp dedicated circuit



Refrigeration specifications

Refrigerant R134a, 11.4 oz (322g)

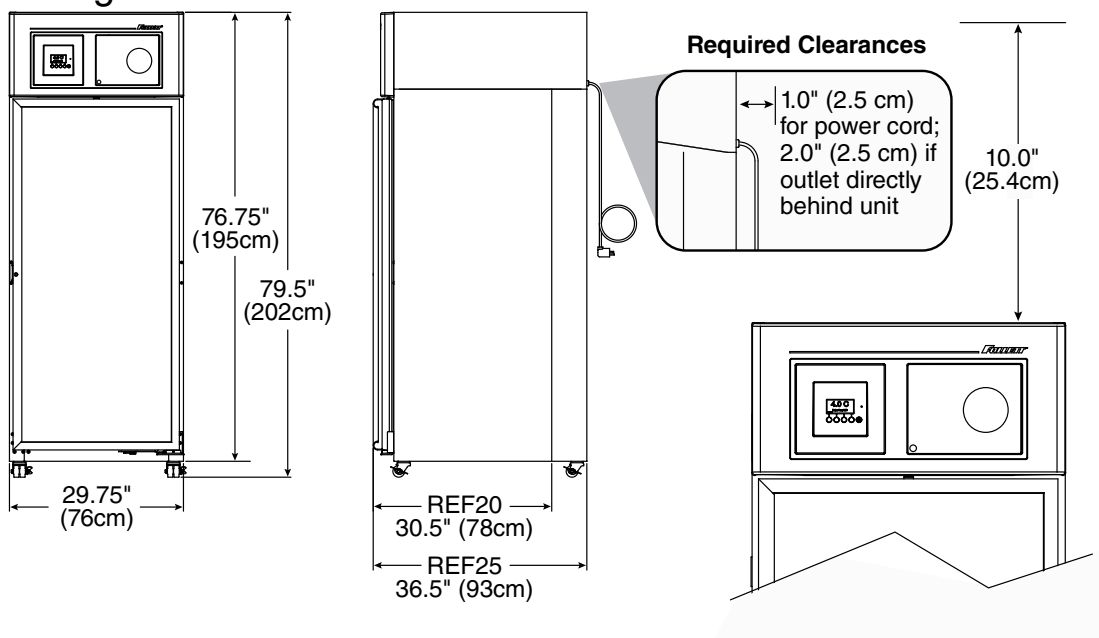Detail drawing




Installation

 DANGER	
	<ul style="list-style-type: none"> • Do not tilt any unit further than 30° off vertical during uncrating or installation • Refrigeration module area contains mechanical, moving parts. Keep hands and arms clear of this area at all times. If access to this area is required, power to unit must be disconnected first.

1. Remove key taped to side of refrigerator cabinet and unlock door.
2. Remove casters from box packed inside refrigerator.
Note: California facilities requiring wall and floor anchors – refer to wall anchor kit instructions (#00938241) packed with seismic kit before proceeding with caster installation.
3. Unbolt refrigerator from pallet.
4. Rotate refrigerator on pallet to install casters (Fig. 1.1).
5. Install casters by hand one at a time using channel locks to tighten, until caster meets against the bottom of cabinet (Fig. 1.2).
6. Peel protective film from exterior walls of stainless cabinet.


Fig. 1

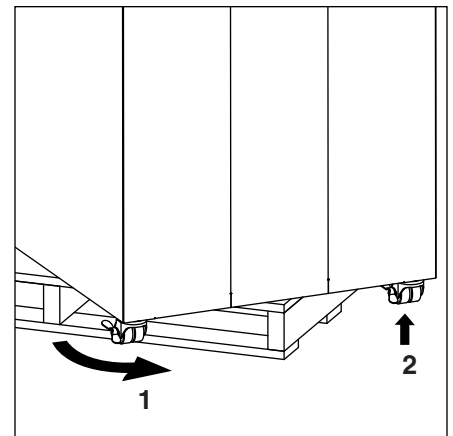

Adjust self-closing door

1. Roll refrigerator to desired location and lock as many casters as possible (Fig 2.1).
2. If the door appears to be sagging, remove the philips screw (Fig. 2.2) holding the Hold Open Bracket to the Hinge Bracket to gain access to the Sag Adjustment Screw (Fig. 2.3). Turn the Sag Adjustment Screw to align the door squarely with the freezer cabinet. When finished, reinstall the Hold Open Bracket.
3. Check door for closing tension and adjust if necessary. To check closing tension, open door 1” and turn screw (Fig. 2.4) counterclockwise until you achieve positive close. (Light will go out) Then, turn another half turn counterclockwise.


Fig. 2

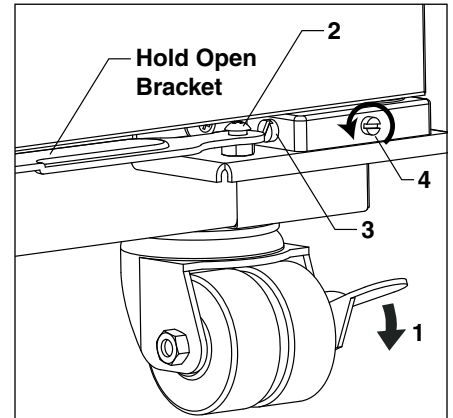

Install shelves (if equipped) and power up

1. If refrigerator is equipped with shelves, remove bag containing shelf supports and position them in desired locations on each pilaster and insert shelves (Fig. 3).
2. Plug refrigerator into a 115V 60Hz 15 amp dedicated outlet.

Fig. 3

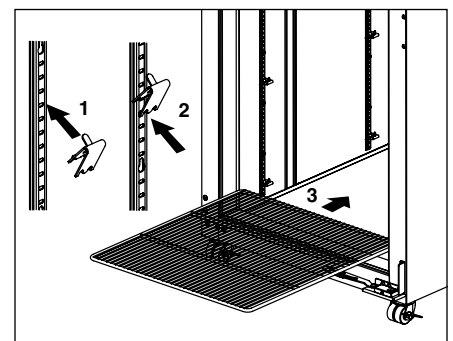

Fill product bottle

1. Remove probe and cap from the top bottle and fill with solution used at your facility (e.g., 50% glycerine and 50% water) (Fig. 4). Replace cap and reinsert probe sensor.
2. If equipped with a bottom probe, perform same procedure for bottom bottle.

ATTENTION INSTALLER
It is imperative that the top probe bottle be filled with a solution that your facility uses to ensure proper operation of unit.

3. Adhere supplied probe wire clips (x4, packaged with manual) to inside wall leading to the probe bottle.
 4. Route probe wire into clips and close the clips.
-
5. Press lower left corner of front facade door to access power switch inside (Fig. 5.1).

6. Press power switch to turn unit on (Fig. 6.1).
7. Install (2) 9V back-up interface batteries (Fig. 6.2) and (1) 9V chart recorder back-up battery (if equipped.)

Fig. 4

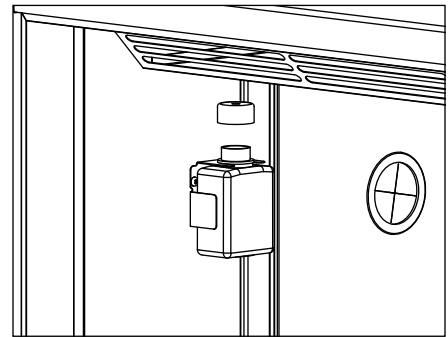


Fig. 5


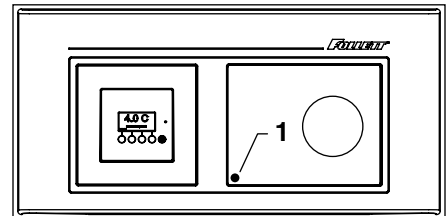


Fig. 6

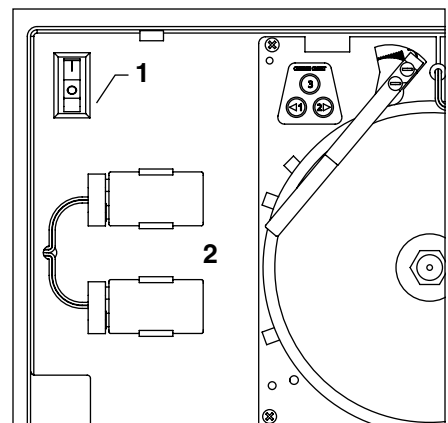

Remote alarm connections (if desired)

Remote communication connections are located on back side of top facade (Fig. 7).


Dry contacts

- Connection to NO or NC contacts provides remote alarming capability
- No temperature data is communicated

RS-485 port

- Connection provides real-time data streaming
 - Temperature
 - Alarms
 - Refrigeration operation
- MODBUS protocol
- Uses 6 pin, 4 conductor RJ11 jack (not supplied)


Fig. 7

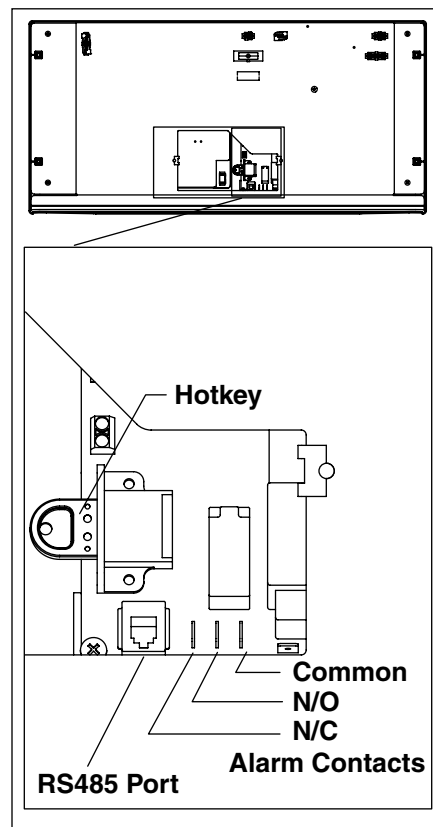

Operation

Quick setup checklist

- Set time/date
- Change temperature display from factory default Celsius to Fahrenheit (if desired)
- Set high/low temperature alarms
- Select password access (if required)
- Setup chart recorder (if equipped)
- Verify solution is in bottle

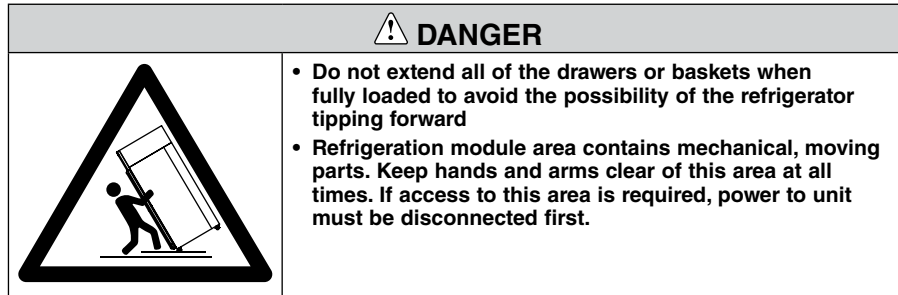


Interior LED lights

Interior lighting of refrigerator with door closed is controlled by light switch (Fig 8.1) on interface.

Lights will be on whenever door is open regardless of light switch selection.

Fig. 8

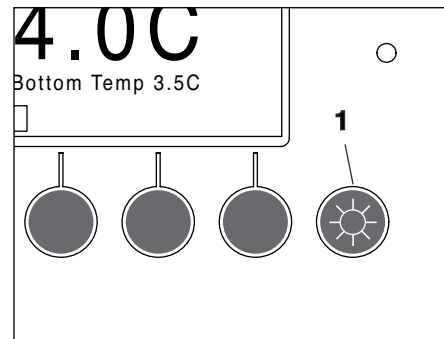

ATTENTION


- The time/date are not retained in the event of a power interruption or if the batteries are dead.
- Change batteries with the unit powered up. Change one battery at a time to retain time/date. Failure to follow this procedure will result in loss of time & date and will not provide accurate time and date stamping for high or low temperature alarms.

Set time & date

Set time and date for stamping of minimum and maximum temperatures.


1. Press **Menu** button (Fig. 9).

Fig. 9

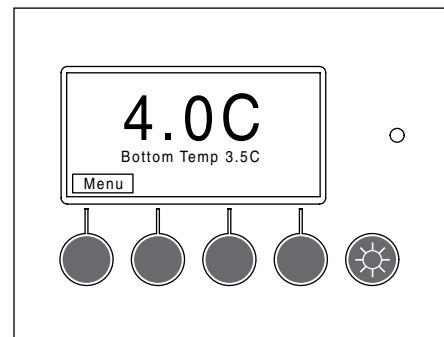


- Use ▲/▼ to scroll and highlight **Set time/date** and press **Enter** to select (Fig. 10).

Fig. 10

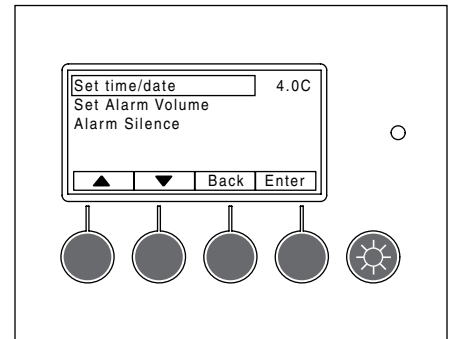

- Adjust the year using ▲/▼ and press **NEXT**.
- Adjust the month using ▲/▼ and press **NEXT**.
- Adjust the day using ▲/▼ and press **NEXT**.
- Adjust the hour using ▲/▼ and press **NEXT**. (24 HR Clock)
- Adjust the minute using ▲/▼.
- Press **Enter** to select values.
- Press **Back** until the main temperature display screen is displayed.

Fig. 11

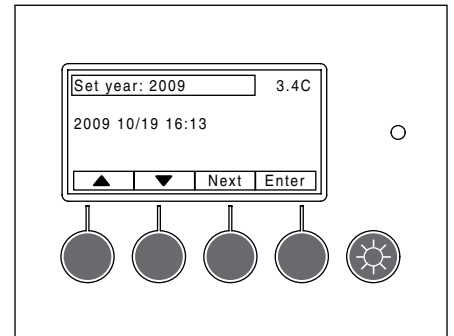

Change temperature display to show degrees Fahrenheit (if desired)

Follett has pre-set the temperature display to read in degrees Celsius. If degrees Fahrenheit display is preferred:

- Press **Menu** button (Fig. 12).
- Use ▲/▼ to scroll and highlight **Display Settings** and press **Enter** to select (Fig. 13).
- Use ▲/▼ to scroll and highlight **Degrees F/C** and press **Enter** to select.
- Use ▲/▼ to scroll and highlight **F** and press **Enter** to select.
- Press **Back** button until the main temperature display screen is displayed. Temperature is displayed in degrees F.

Fig. 12


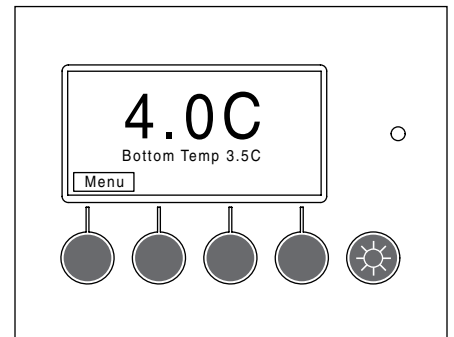


Fig. 13

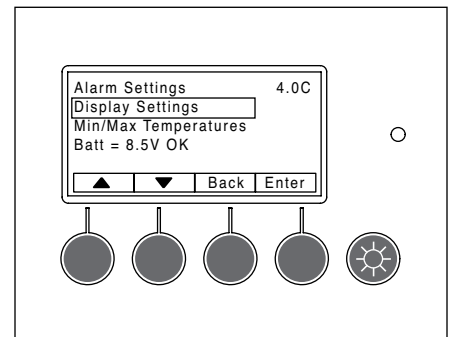


High and low alarm setup

The high alarm is set to a default of 40 C (104 F) from the factory. This default setting is used so that the alarm will not sound while the unit is pulling down to temperature. The low alarm default setting is 1.0 C (33.8 F).

After the unit has been installed and the system has stabilized (approximately 1 hr), the high and low alarm set points should be adjusted to desired settings.


1. Press **Menu** button (Fig. 14).

Fig. 14

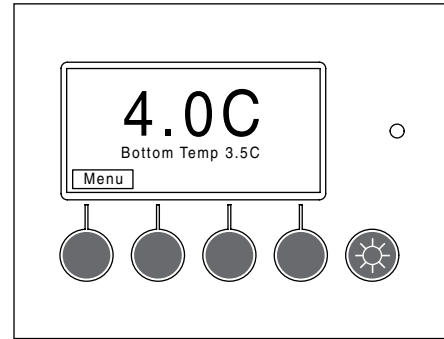


2. **Alarm Settings** will be highlighted. Press **Enter** to select (Fig. 15).

Fig. 15

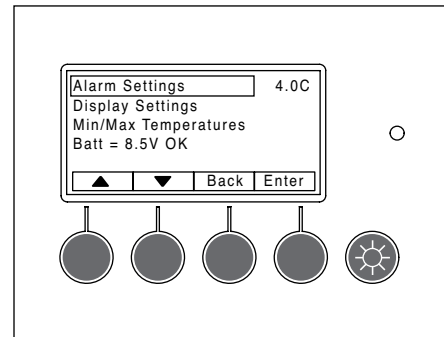


3. Use ▲/▼ to scroll and highlight **High Alarm Set** and press **Enter** to select (Fig. 16).

Fig. 16

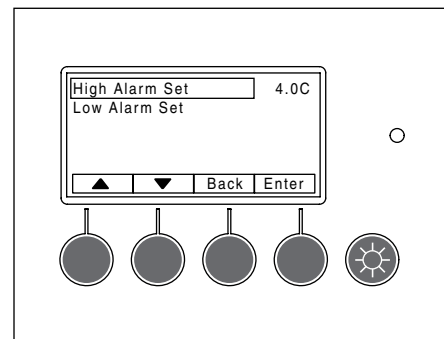


4. Use ▲/▼ to change set point and press Enter to accept (Fig. 17).

Fig. 17

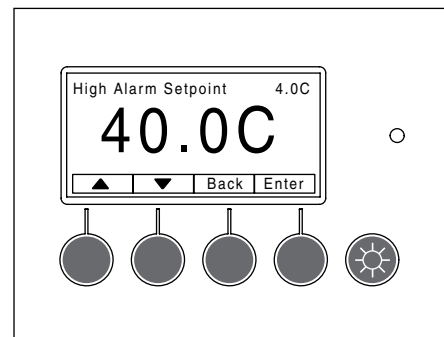


5. Use ▲/▼ to scroll and highlight **Low Alarm Set** and press **Enter** to select (Fig. 18).

Fig. 18

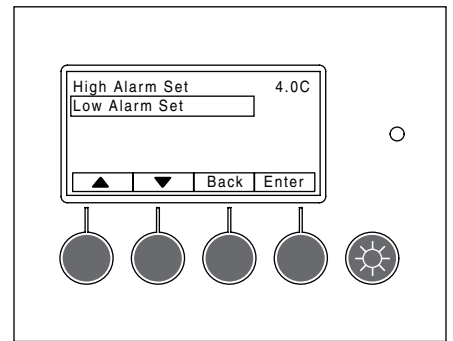

6. Use ▲/▼ to change set point and press Enter to accept (Fig. 19).
7. Press **Back** until the main temperature display screen is displayed.

Fig. 19

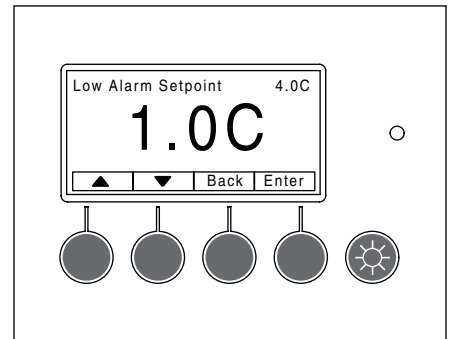

Alarms

The refrigerator has several alarms that will sound an audible and visual alarm. Alarm type is shown on the top line of the alarm display and multiple alarms will show progressively on the display. A red LED, to the right of the display, accompanies all alarms.


All active alarms are conveyed to remote alarm relays. Remote alarm relay provides a choice of either a normally closed (NC) or normally open (NO) contact. It is recommended for use at 24V 1A (relay is rated for 250V 10A).

Alarm displayed	Conditions
Temp Alarm	Only top product temperature probe activates this alarm. Bottom product temperature probe, if installed, is for display reference only.
Top Probe Error	Activates with loss of signal from top probe (open or short circuit).
Bottom Probe Error	Activates with loss of signal from bottom probe (open or short circuit).
Door Ajar Alarm	Activates with door open longer than 1 minute.
AC Power Alarm	Activates after 2 min delay with loss of AC power. Display backlight will turn off to conserve battery life, but pressing any button will brighten display for 30 seconds.
Low Battery Alarms	Display activates when (2) 9V controller back-up batteries decrease to 7.5V (maximum 4 hr back-up time remaining.) This alarm will audibly chirp every 5 minutes. Chart recorder – LED flashes green. All batteries should be changed at same time. Replace batteries one at a time with AC power on to keep correct time/date.
Evap Probe Error	Activates with loss of signal from refrigeration control evaporator probe (open or short circuit). Refrigerator goes into a fixed on/off cycle until corrected. Default settings are 4 minutes compressor on time and 8 minutes compressor off time. Cycle on time and off time are adjustable in faulty probe cycle sub-menu of refrigeration controls menu.

Alarm muting

All audible alarms can be muted for up to 60 minutes in 5 minute increments by pressing the **MUTE** button. The desired mute period will display, and count down, in the window adjacent the MUTE label (Fig. 20).

Fig. 20

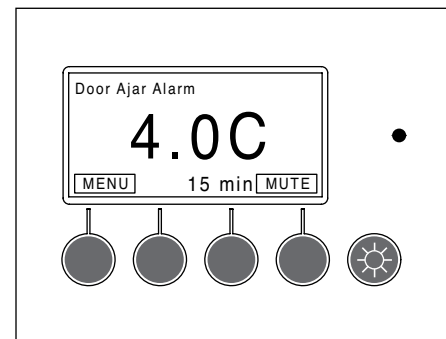


Alarm volume selection

Factory default setting is “high”. To change volume of the alarm:


1. Press **Menu** button (Fig. 21).

Fig. 21

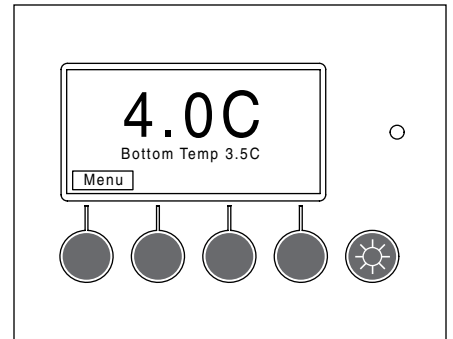


2. Use ▲/▼ to scroll and highlight **Set Alarm Volume** and press **Enter** to select (Fig. 22).

Fig. 22

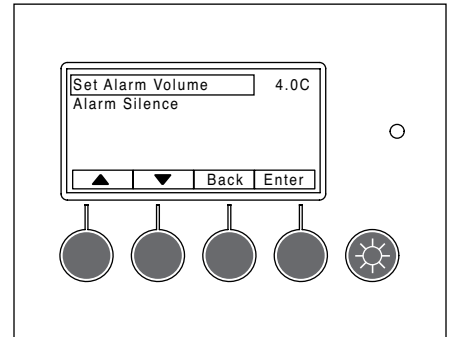

3. Use ▲/▼ to select the volume (high/medium/low) and press **Enter** to select (Fig. 23).
4. Press **Back** until the main temperature display screen is displayed.

Fig. 23

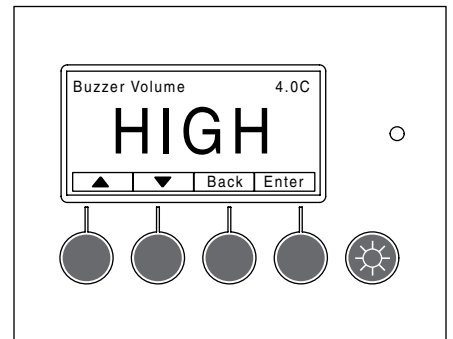


Alarm silence

Audible alarming can be turned off through the menu interface:


1. Press **Menu** button (Fig. 24).

Fig. 24

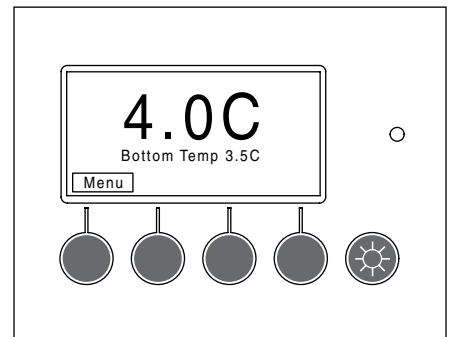

- Use ▲/▼ to scroll and highlight **Alarm Silence** and press **Enter** to select (Fig. 25).

Fig. 25

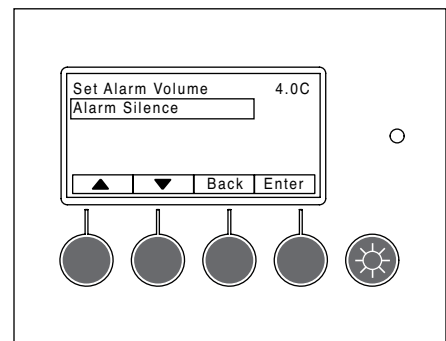

- Use ▲/▼ to select **ON** or **OFF** and press **Enter** to select (Fig. 26).
- Press **Back** until main temperature display screen is displayed.

Fig. 26

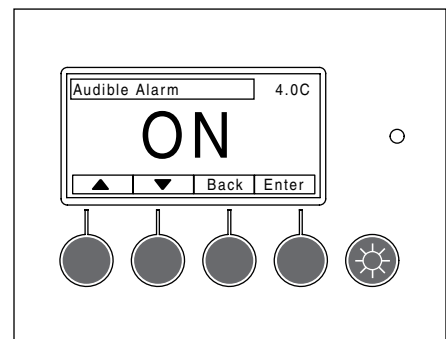


Fig. 27


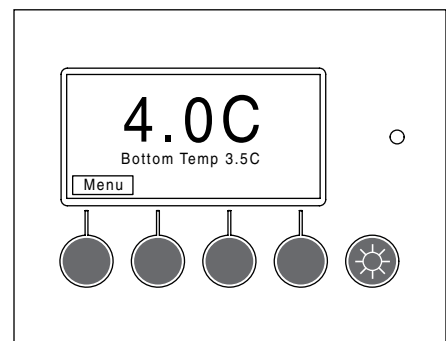


Fig. 28

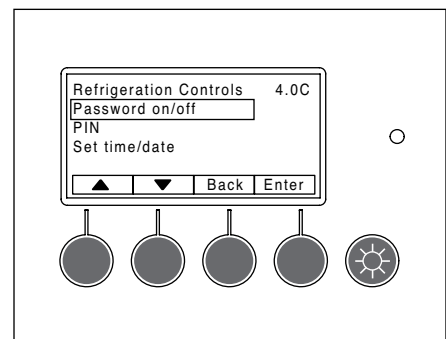

Password PIN security (default is 3843)

If desired, access to the user selectable features and settings can be controlled by a password PIN. The refrigerator ships with the password option turned off. The default password PIN, if the password is turned on, is 3843. This password pin number will always work.

To turn password on/off

- Press **Menu** button (Fig. 27).
- Use ▲/▼ to scroll and highlight **Password on/off** option and press **Enter** to select (Fig. 28).

3. Use ▲/▼ to choose **ON** or **OFF** and press **Enter** to select (Fig. 29).
4. Press **Back** until the main temperature display screen is displayed.

Fig. 29

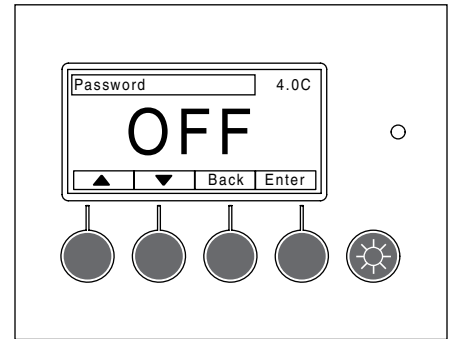


Fig. 30

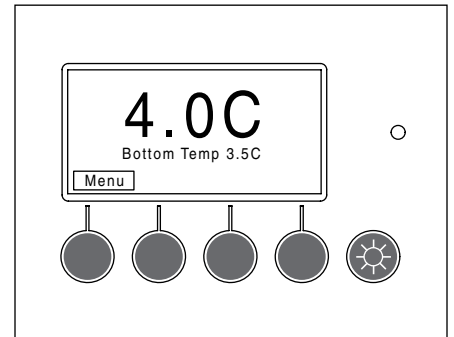


Fig. 31


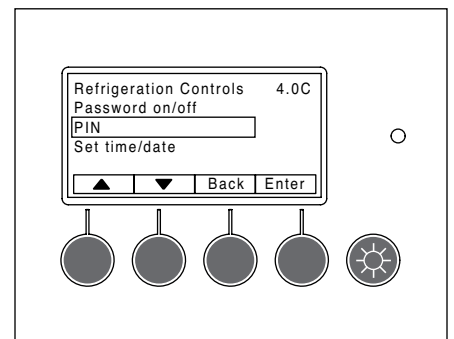


Fig. 32

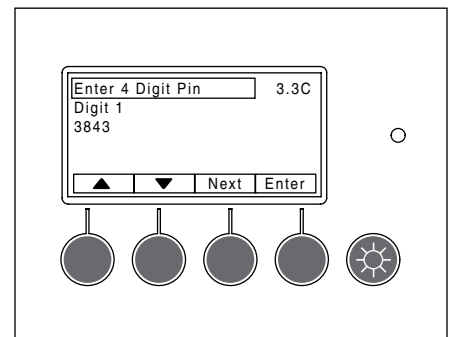

To make a new password PIN


1. Press **Menu** button (Fig. 30).
2. Use ▲/▼ to scroll and highlight **PIN** and press **Enter** to select (Fig. 31).
3. Use ▲/▼ to select the first number of the 4 digit pin. Use the **Next** button to move to the next digit, adjust with ▲/▼ and repeat for the last 2 digits. Press **Enter** to select (Fig. 32).
4. Press **Back** until the main temperature display screen is displayed.

Min/max temperature logging

The controller retains maximum and minimum temperatures since last reset and provides a date and time stamp of when they occurred.


1. Press **Menu** button (Fig. 33).

Fig. 33

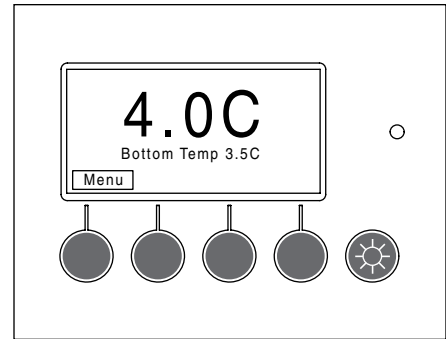


2. Use ▲/▼ to scroll and highlight **Min/Max Temperatures** and press **Enter** to select (Fig. 34).

Fig. 34

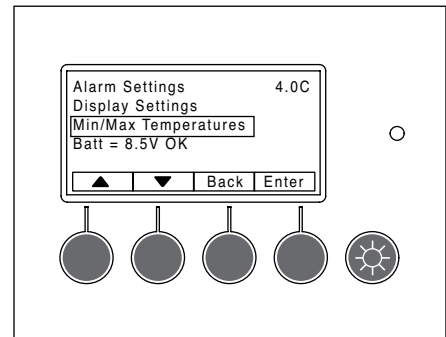

3. Use ▲/▼ to select the min or max logged temperature (Fig. 35). Press **Enter** to reset log.
4. Press **Back** until the main temperature display screen is displayed.

Fig. 35

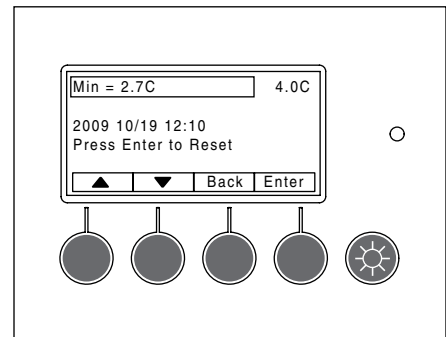


Probe calibration

Follett refrigerator probes are calibrated to a known standard prior to leaving the factory. Recalibration can be easily accomplished at any time, if desired:


1. Remove upper temperature probe from probe bottle and completely submerge into an ice bath.
Note: Ice bath should be densely compacted ice (flake, crushed, nugget, or cracked) and a little water.
2. Place ice bath with probe in refrigerator and close door.
3. Press **Menu** button (Fig. 36).

Fig. 36

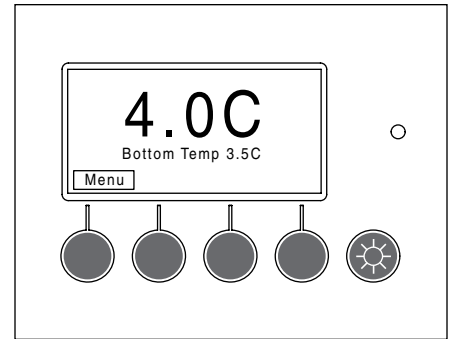


4. Use ▲/▼ to scroll and highlight **Refrigeration Controls**, press and hold **Enter** and ▼ together for 3 seconds (Fig. 37).

Fig. 37

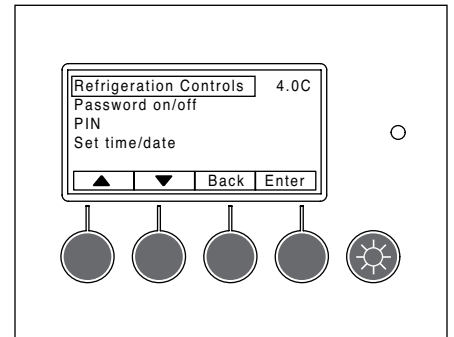


5. **Calibration** should be highlighted, press **Enter** to select (Fig. 38).

Fig. 38

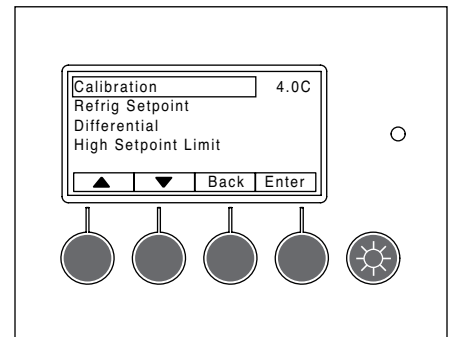

6. **Cal Top Probe** should be highlighted, press **Enter** to select (Fig. 39).

Fig. 39

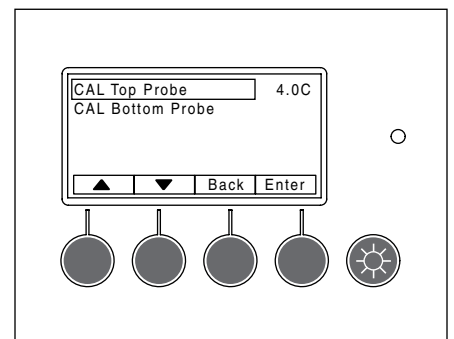

7. When probe temp stabilizes press **Enter** button calibrating it to 0.0 C (32 F) (Fig. 40). Wait one minute and press **Enter** button again to confirm calibration.
8. Remove probe from ice bath and return to solution-filled probe bottle.
9. Press **Back** until main temperature display screen is displayed.
10. Repeat steps for bottom probe, if equipped.

ATTENTION
<ul style="list-style-type: none"> • Bottom temperature probe is only standard on blood bank models • If unit has the chart recorder option, refer to “Chart recorder calibration” on page 18.

Fig. 40

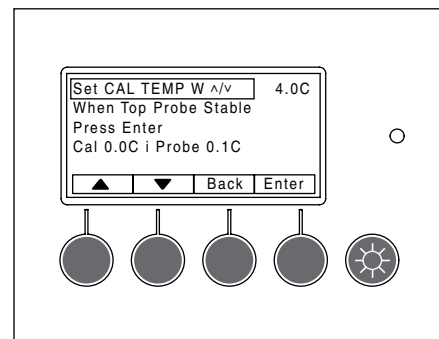


Chart recorder set up (if so equipped)

The chart recorder is mounted in the front right of the facade behind the door (Fig. 41). A package of 50 charts that record in C is included with your refrigerator. If you have changed your display to show temperatures in F, you will need different charts. Please call Follett at (877) 612-5086 or +1 (610) 523-9361 to order part number 00967729.

The chart recorder is powered by the controller and has its own battery back up. An LED provides battery status info: Steady green = battery okay, Flashing green = low battery or no power to the controller.

Chart recorder calibration

If desired, calibration should be done at the same time as probe calibration (page 17) by pressing either the left or right arrow keys to correspond with the calibrated product display temperature.

1. Press **Left #1** or **Right #2** arrow button (Fig. 41) for 5 seconds until pen begins to move.
2. Continue to press **#1** or **#2** button to move pen to record same temperature as probe.


Changing the chart paper

1. Press and hold **#3** button until pen begins to move off chart.
2. Once the pen moves completely off chart, unscrew knob (counter-clockwise) at center of chart.
3. Remove old chart and place new chart in position so that correct time coincides with time line groove on recorder.
4. Reinstall knob.
5. Press and hold **#3** button for approximately (1) one second until pen begins to move back onto chart.
6. Make sure pen is marking paper. If not, carefully adjust arm so that pen makes contact with paper.

Replacing the pen

1. To replace pen, press and hold the **#3** button until pen starts to move off chart.
2. Once pen has moved completely off chart, carefully lift pen up and unsnap the “U” clip part to remove pen from metal pen arm.
3. Position new pen on metal pen arm and snap “U” clip underneath pen around arm to secure.
4. Press and hold **#3** button for approximately (1) second until pen begins to move back onto chart.
5. Make sure pen is marking. If not, carefully adjust arm so that pen makes contact with paper.


Fig. 41

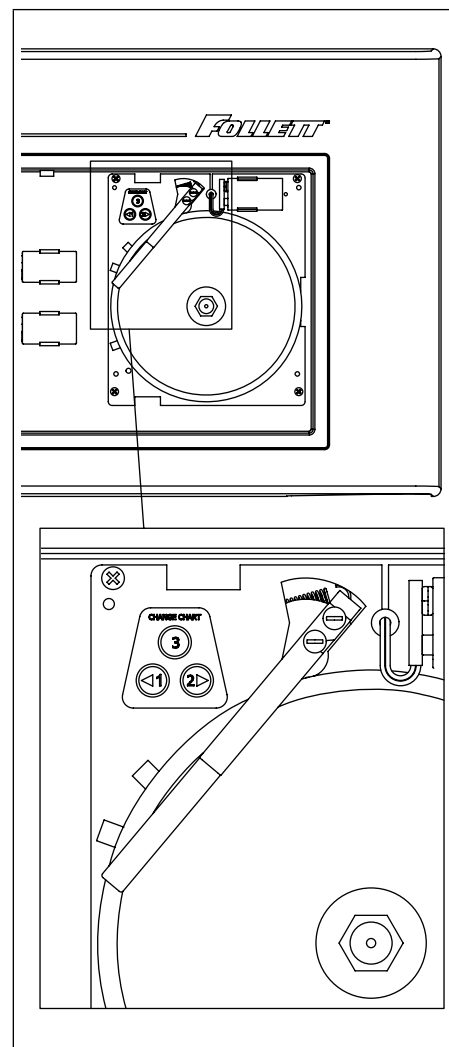

Removing drawers

1. Pull drawer forward to stop (Fig. 42.1).
2. Lift drawer front to free front rollers from slides (Fig. 42.2).
3. Still lifted, pull drawer forward to free back rollers from slides.

Fig. 42

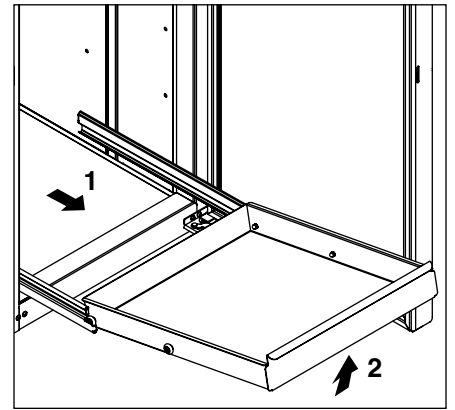

Removing slides

1. Push slides all the way back.
2. Swing bottom of slide away from refrigerator wall and lift slide off rollers.

Cleaning

Use non-chlorine-based cleaners. Cleaners containing chlorine can cause staining and pitting of the stainless steel.

Disconnect power to unit by turning the power switch off, located on the upper front facade panel, and removing the power cord from the receptacle.

Interior cleaning

Using a sponge or soft cloth, clean unit with a non-abrasive, non-chlorinated, all-purpose detergent. Note: The left and right air plenums should be removed to clean them, as well as behind them (Fig. 43, 44). In addition the shelves, baskets, drawers and slides should be cleaned as well.

Exterior cleaning

Wipe stainless steel exterior with a soft cloth in the direction of grain as needed. Stainless steel polish may be used to enhance the finish of the unit. The glass door and exterior parts may be cleaned with a soft cloth, window cleaner or other non-abrasive cleaner.

Fig. 43


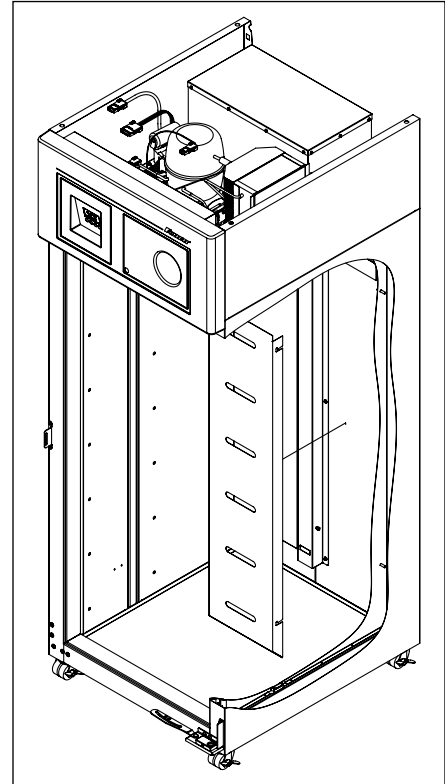


Fig. 44

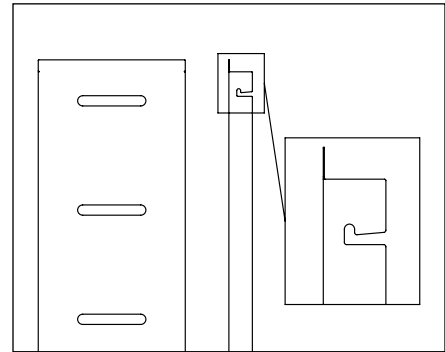


Annual cleaning

Removal of dust and other particulates from air intake areas and the condenser is important for proper operation. Environments with large amounts of dust may require more frequent cleaning.


1. Disconnect power to unit by turning power switch off (located on upper front facade panel) and removing power cord from receptacle.
2. The condenser can be accessed from top for cleaning. The side panel in front of condenser may be completely removed for easier cleaning by disconnecting facade and removing upper side panel. See “Facade removal” on page 27 for facade removal instructions.
3. Inspect drain pan for any debris or obstruction in condensate pan (Fig. 45.1).
4. Use a vacuum cleaner with brush attachment to clean condenser, compressor and related parts.
5. Reinstall upper side panel and facade if removed.

Fig. 45

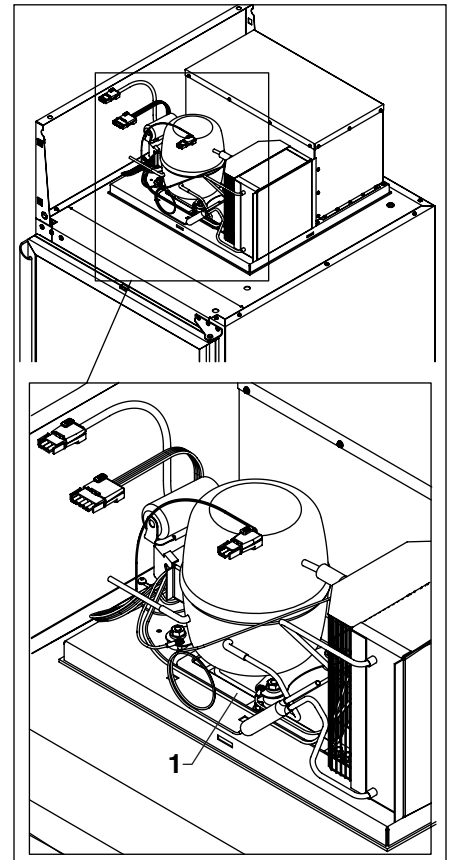

6. Door maintenance: Place a dab of white lithium grease on top of hold open arm and keeper on lower hinge assembly of door (Fig. 46.1).

Fig. 46

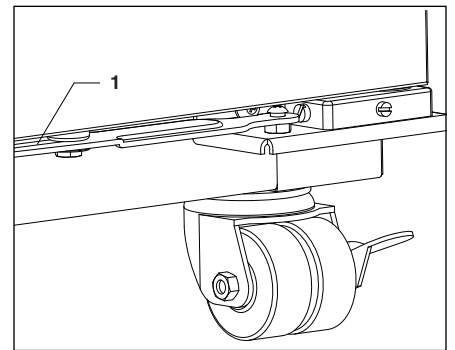

Service

The temperature controller and evaporator probe indicate when the refrigeration system is required to turn on and off.

The refrigeration system removes heat from the cabinet interior and rejects it to the surrounding room air. When the evaporator probe temperature reaches 5.0 C (9.0 F) above the controller set point, the probe signals the controller to turn the refrigeration system on. The normally open controller contacts close and energize the condenser fan motor and compressor. The compressor uses a current-style starting relay and a starting capacitor to start the compressor motor.


When the evaporator temperature falls to 5.0 C (9.0 F) below the set point, the probe signals the controller to turn the refrigeration system off. The controller contacts reopen, which de-energizes the condenser fan motor and the compressor. The evaporator fan motor runs continuously.

Any accumulated frost on the evaporator coils melts during the off cycle. The condensate drains to a drain pan mounted in the base of the refrigeration unit. The pan is located under the compressor. The heat from the compressor, condenser and fan evaporates any condensate in the drain pan as well as providing preliminary cooling of the high pressure vapor prior to entering the condenser.

Refrigeration system

The REF20/25 series refrigeration system is designed to give many years of trouble-free service. Except for routine cleaning of the air-cooled condenser and related parts, the refrigeration system requires no service or maintenance. The system uses a capillary tube metering device and is critically charged. Access fittings are not provided. Temporary piercing valves are required to work on the refrigeration system. Follett recommends that if hoses are ever connected to the refrigeration system for service, the refrigerant should be recovered, permanent access valves should be installed and the system evacuated, and recharged by weighing in the correct refrigerant charge (refrigerant R134a, 11.4 oz (322g).

Note: Do not charge the system by pressures.

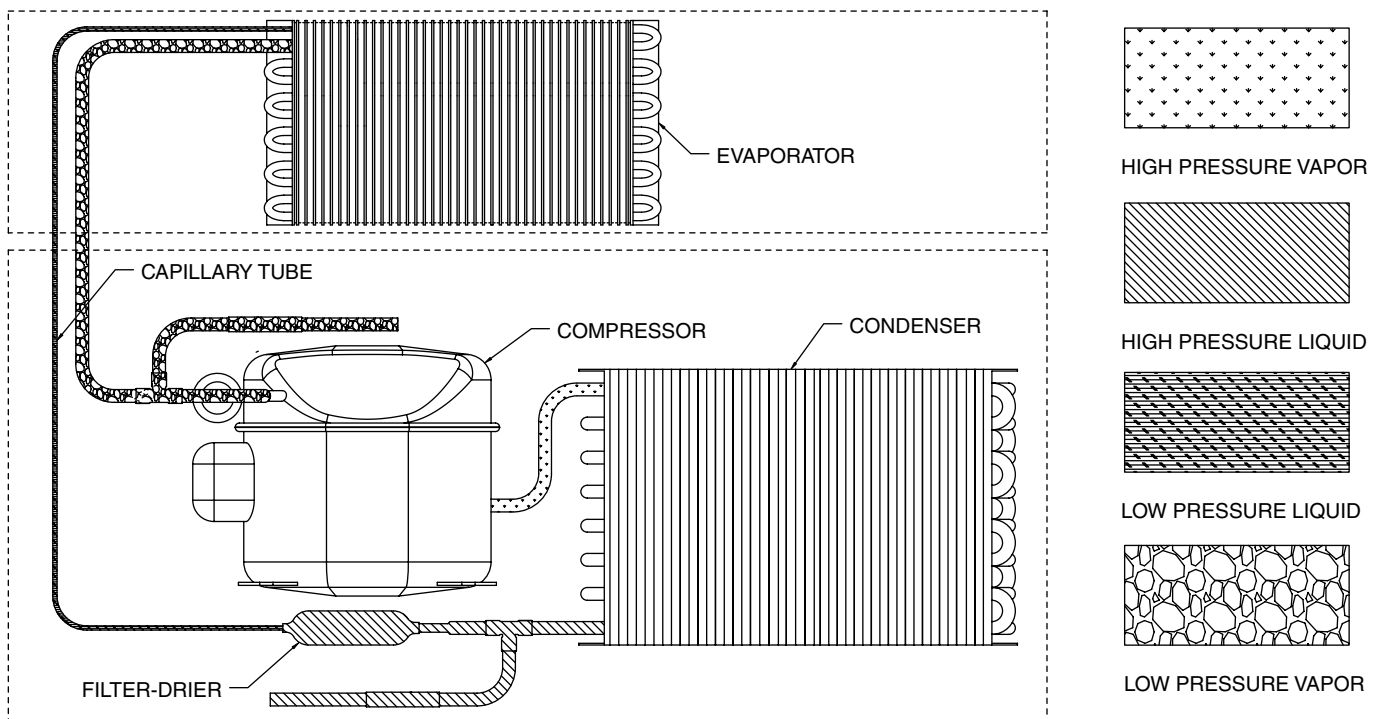

Defrosting

REF20/25 series upright refrigerators do not require manual defrosting. The evaporator defrosts automatically when the refrigeration unit is in the OFF cycle. The evaporator fan will run continuously.

Controller operation

ATTENTION
BEFORE MAKING ANY CHANGES TO FACTORY-DEFAULT REFRIGERATION SETTINGS, CONTACT FACTORY.

The controller displays product temperatures in degrees C or F (C is factory default) to 1 decimal point.

The controller is pre-programmed with a 0.5 C (32.9 F) set point which provides a compressor cut-in at 5.5 C (41.9 F) evaporator coil temperature and cut-out at -4.5 C (23.9 F) evaporator coil temperature. This will correlate to a product temperature of approximately 4.0 C (39.2 F).

All set points have a 5 C (9.0 F) differential. The 5.0 C (9.0 F) differential means that with a 0.5 C (32.9 F) set point, for example, the compressor will turn off at -4.5 C (23.9 F) evaporator temperature and turn on when the evaporator temperature reaches 5.5 C (41.9 F).


Product temperature will be displayed on the digital temperature display and chart recorder (if equipped).

Temperature sensor readings

To view temp sensor probe readings:


1. Press **Menu** button (Fig. 47).

Fig. 47

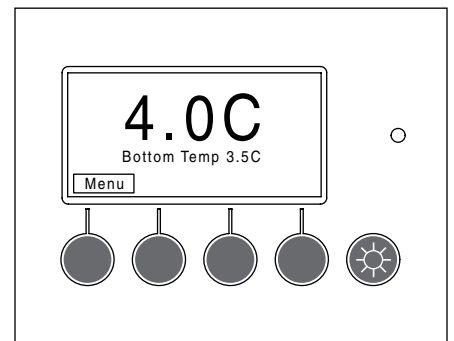


2. Use ▲/▼ to scroll and highlight **Refrigeration Controls** (Fig. 48).

Fig. 48

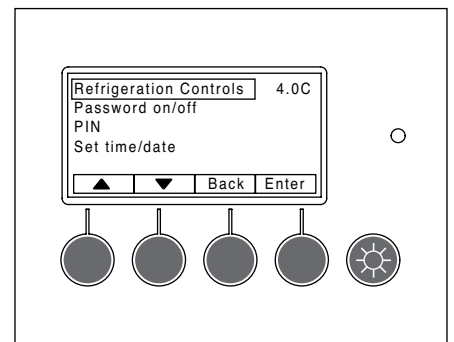


3. Press and hold **Enter** and ▼ together for 3 seconds (Fig. 49).

Fig. 49

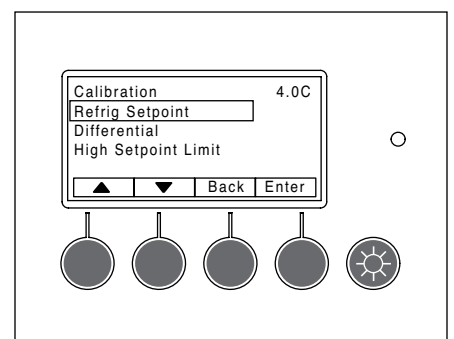


- Use ▲/▼ to scroll and highlight **Temp sensors** and press **Enter** to select (Fig. 50).

Fig. 50

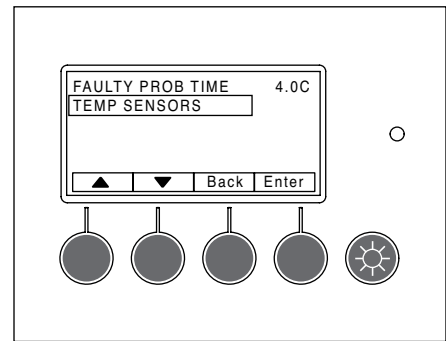

- The following will be displayed (Fig. 51).
 RTD1= Top probe sensor reading
 RTD2= Bottom probe sensor reading
 (if equipped, otherwise displays default 25 C (77 F))
 NTC1= Evaporator Probe reading
 NTC2= Displays a default value and is not used
- Press **Back** until main temperature display screen is displayed.

Fig. 51

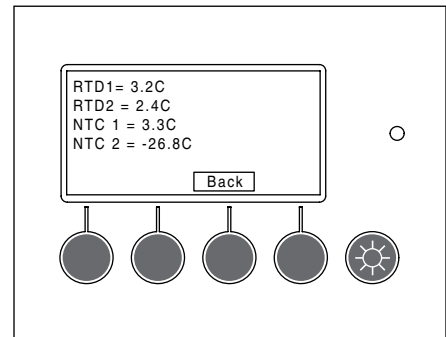


Heated door (option) – changing cycle

Factory default for the heated door and fascia heater cycle is set to 50%, based on a 20 minute cycle (e.g. 50% = 10 minutes off, 10 minutes on).


- Press **Menu** button (Fig. 52).

Fig. 52

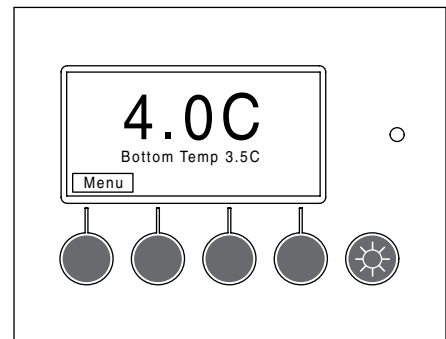


- Use ▲/▼ to scroll and highlight **Refrigeration Controls** (Fig. 53).

Fig. 53

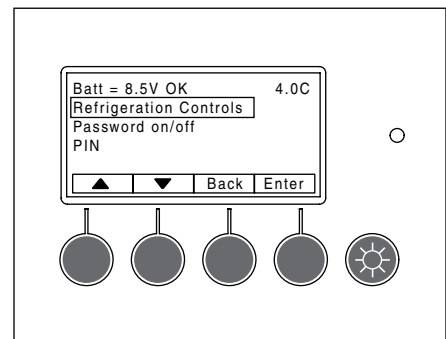


3. Press and hold **Enter** and ▼ together for 3 seconds (Fig. 54).

Fig. 54

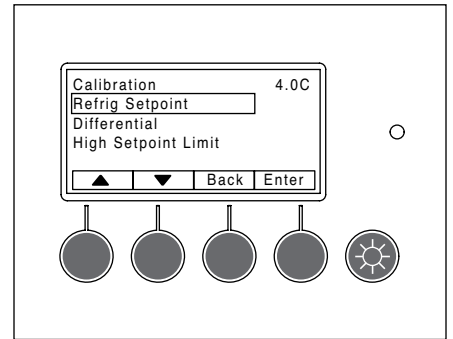


4. Use ▲/▼ to scroll and highlight **Door Heater Cycle** and press **Enter** to select (Fig. 55).

Fig. 55

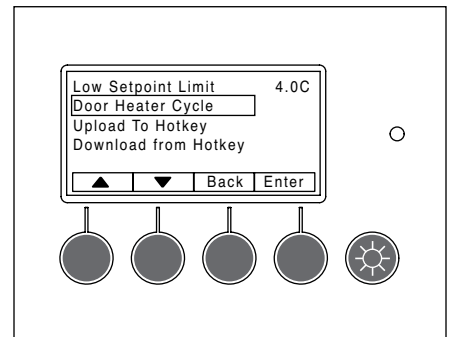

5. Heater ON cycle is shown as a percentage of 10 minute cycle (Fig. 56).
6. Use ▲/▼ to adjust heater ON cycle percentage and press **Enter** to select.
7. Press **Back** until the main temperature display screen is displayed.

Fig. 56

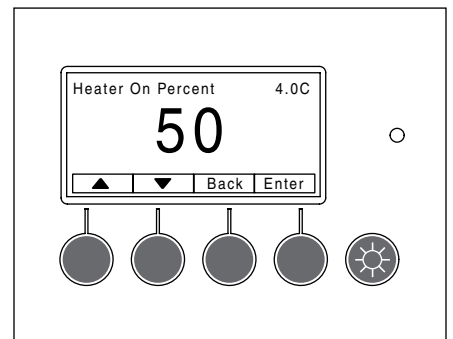

Changing temperature set point

ATTENTION

Before making any changes to factory-default refrigeration settings, contact factory. Factory default is in degrees C. If user has changed display to show temperature in degrees f, all other temperatures will display in degrees F.

The temperature control system is preset by the factory to provide a product temperature of approximately 4.0 C (39.2 F). If desired, the evaporator set point temperature can be changed to achieve a different product temperature. The factory preset 5.0 C (9 F) differential will be maintained regardless of the controller set point.

EXAMPLE:

If the product temperature is 4.0 C (39.2 F) and the product temperature needs to be 3.0 C (37.3 F). Adjust the refrigeration set point down 1.0 C to -0.5 C (1.8 F to 31.1 F) from default setting of 0.5 C (32.9 F).

Changing the temperature set point:

ATTENTION

Cabinet temperature adjustments can be made by simply modifying the temperature control setpoint. Modifications to the factory default settings for the control differential or evaporator offset parameters will result in excessive compressor cycles that may lead to premature compressor/start component failures. Component failures caused by adjusting these critical control operating parameters may void the equipment warranty.

1. Press **Menu** button (Fig. 57).
2. Use ▲/▼ to scroll and highlight **Refrigeration Controls** (Fig. 58).
3. Press and hold **Enter** and then ▼ together for 3 seconds.
4. Scroll ▼ to **Refrig Setpoint**, press **Enter** to select (Fig. 59).

Fig. 57

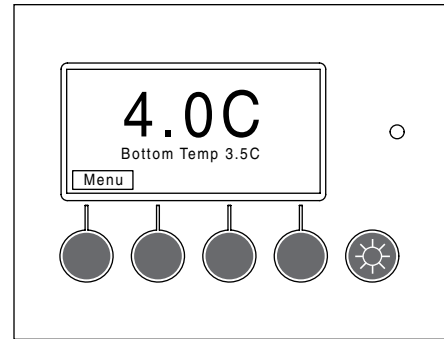


Fig. 58



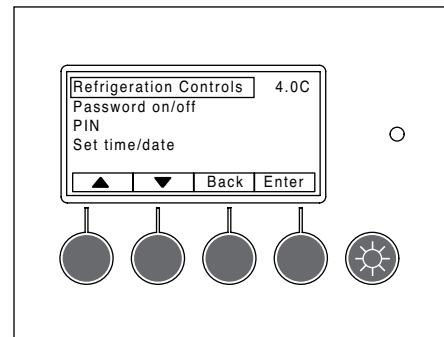


Fig. 59

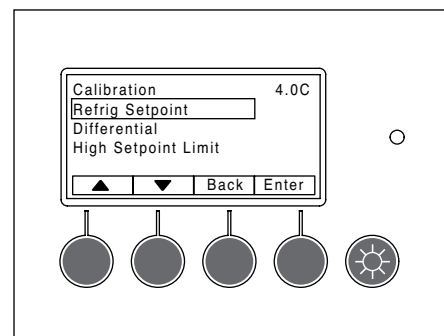

5. Use ▲/▼ to increase or decrease the set point and press **Enter** to select (Fig. 60).
6. Press **Back** until the main temperature display screen is displayed.


Fig. 60

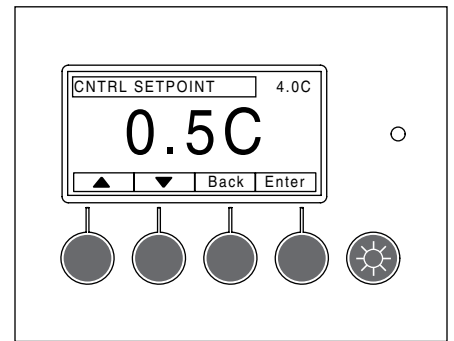



Facade removal

The facade is removed by loosening the top left and right Phillips screws on the back of the facade to allow the securing clips to rotate away from the side panels. The facade can simply pull off of the front by pulling it out of the bayonet mounts that secure it to the upper left and right panels. If the facade will be completely removed, the electrical plugs should be disconnected. The ground wire should be removed from the top right grounding screw on the cabinet box, not from the facade (Fig. 61.1).

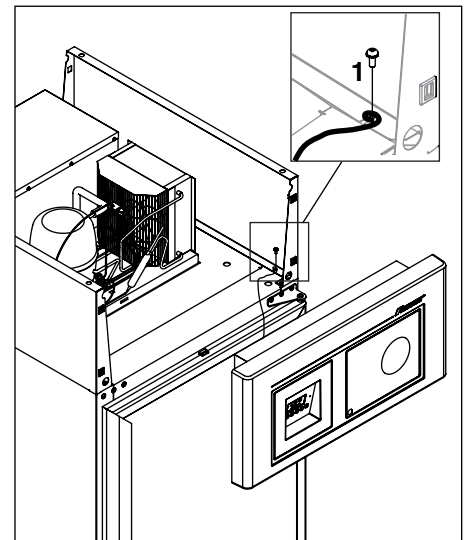Fig. 61




Controller replacement

 CAUTION	
	<ul style="list-style-type: none">• Controller board is susceptible to electrostatic discharge. Extreme care should be exercised by using a grounding strap when handling and installing control board.

1. Turn off power to unit and unplug.
2. Remove facade and disconnect electrical plugs.
3. Remove screw securing ground screw to refrigerator cabinet (Fig. 62.1). Do not remove ground wire from facade.
4. Place facade face down on flat clean surface with bottom of facade facing you, and remove four screws attaching back panel to facade.
5. Lift and rotate back panel toward you.
6. Take precautions for proper grounding to reduce risk of electrostatic discharge (ESD) to the controller board.
7. Remove 5 screws securing board to facade (Fig. 62.2).
8. Reposition board away from mounting location and install new board.
9. Using wiring schematic (attached to inside of removed back panel) as a guide, remove wires one at a time from back of existing controller and connect to corresponding terminals on replacement controller.
10. If ribbon cable from the display is not connected, carefully connect to corresponding location on board (Fig. 62.3).
11. Make sure wiring is secured at same factory locations, so that cover will not pinch any wires when reinstalled.
12. Flip facade cover back into position on facade and reinstall 4 screws.
13. Place facade on top of refrigerator above compressor.
14. Reconnect electrical plugs and reinstall screw for grounding wire to grounding location on cabinet.
15. Remount facade to bayonet mounts.

Fig. 62

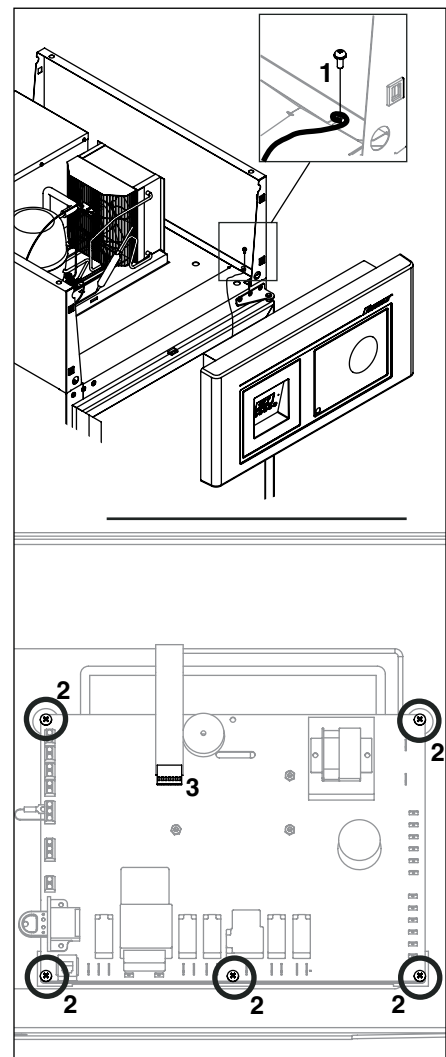

Reversing the door swing (requires a new door)

REF20/25 units ship standard with a non heated right hinged door. To change to a left hinged configuration requires a new door. For heated door options other parts may be required. If the unit has a heated door option consult technical service by calling (877) 612-5086 or +1 (610) 252-7301.


Heated doors

Heated doors are factory installed. An unheated door cannot be retrofitted to a heated door in the field.

Door gasket replacement

1. Remove existing gasket from mounting track (Fig. 63).
2. Verify mounting track is free of any remaining gasket material.
3. Align new gasket with mounting track and press firmly in place.
4. Open and close door, checking for proper gasket seal without pinching against refrigerator.

Fig. 63

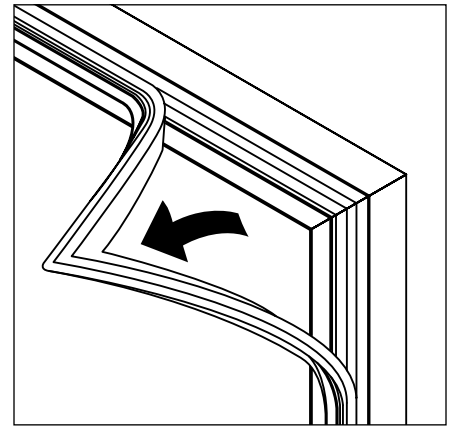


Slide-out unit cooler


Follett's slide-out refrigeration system allows technicians to remove the entire unit from the refrigerator top.

1. If necessary remove facade and upper side panels.
2. Remove four bolts securing unit cooler to refrigerator top (Fig. 64).
3. Gently lift off cooler.


Fig. 64

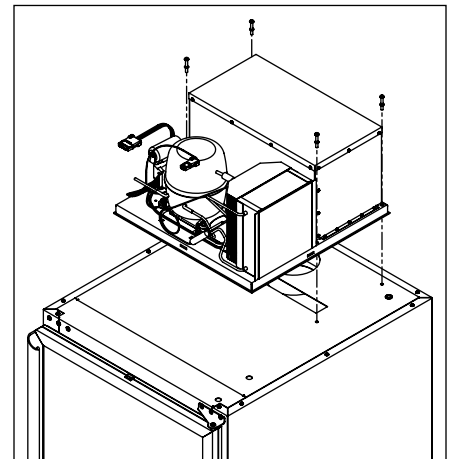

Wiring diagram (After Serial Number E15526)

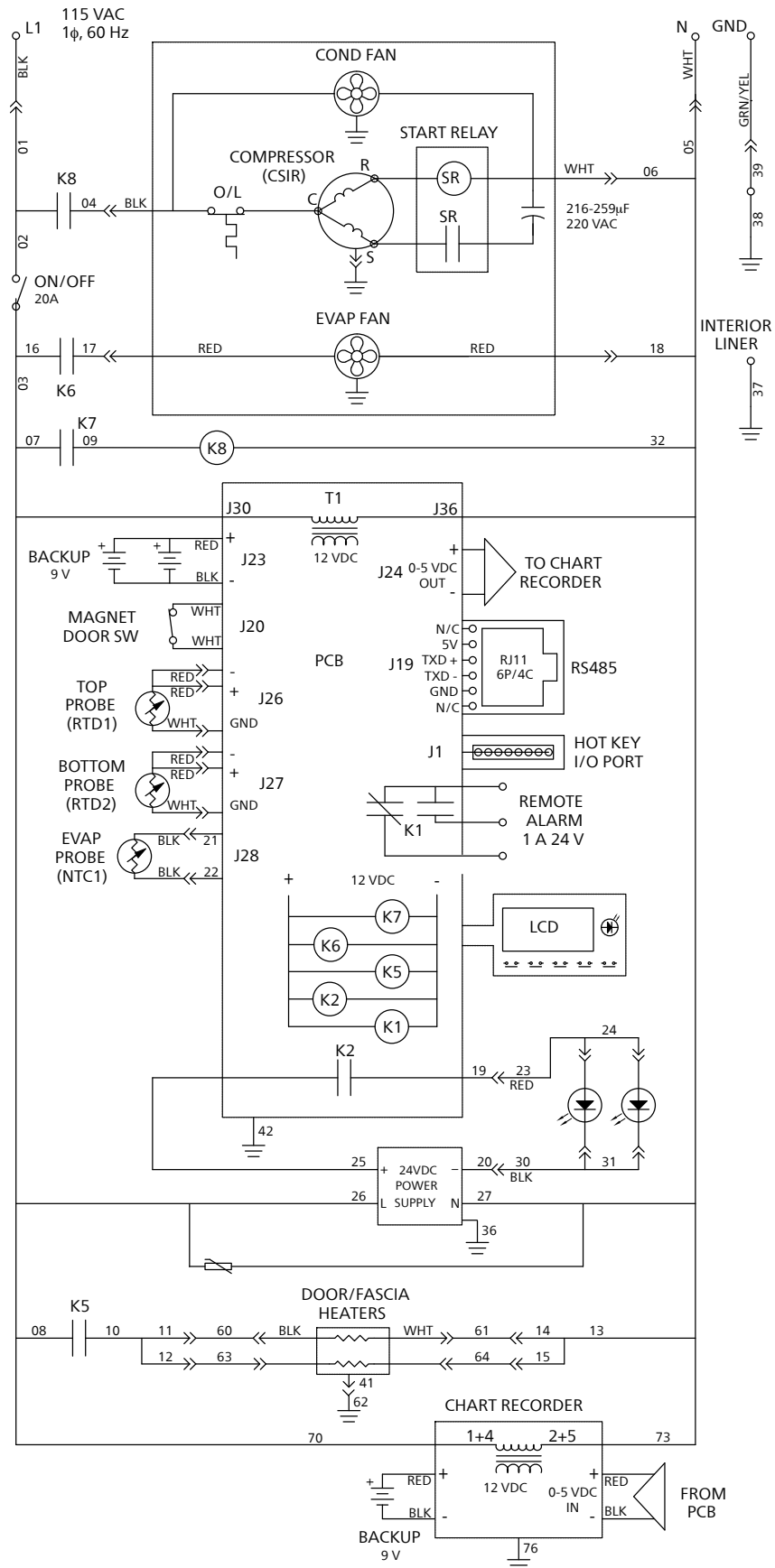

Wiring diagram (After Serial Number E15526)

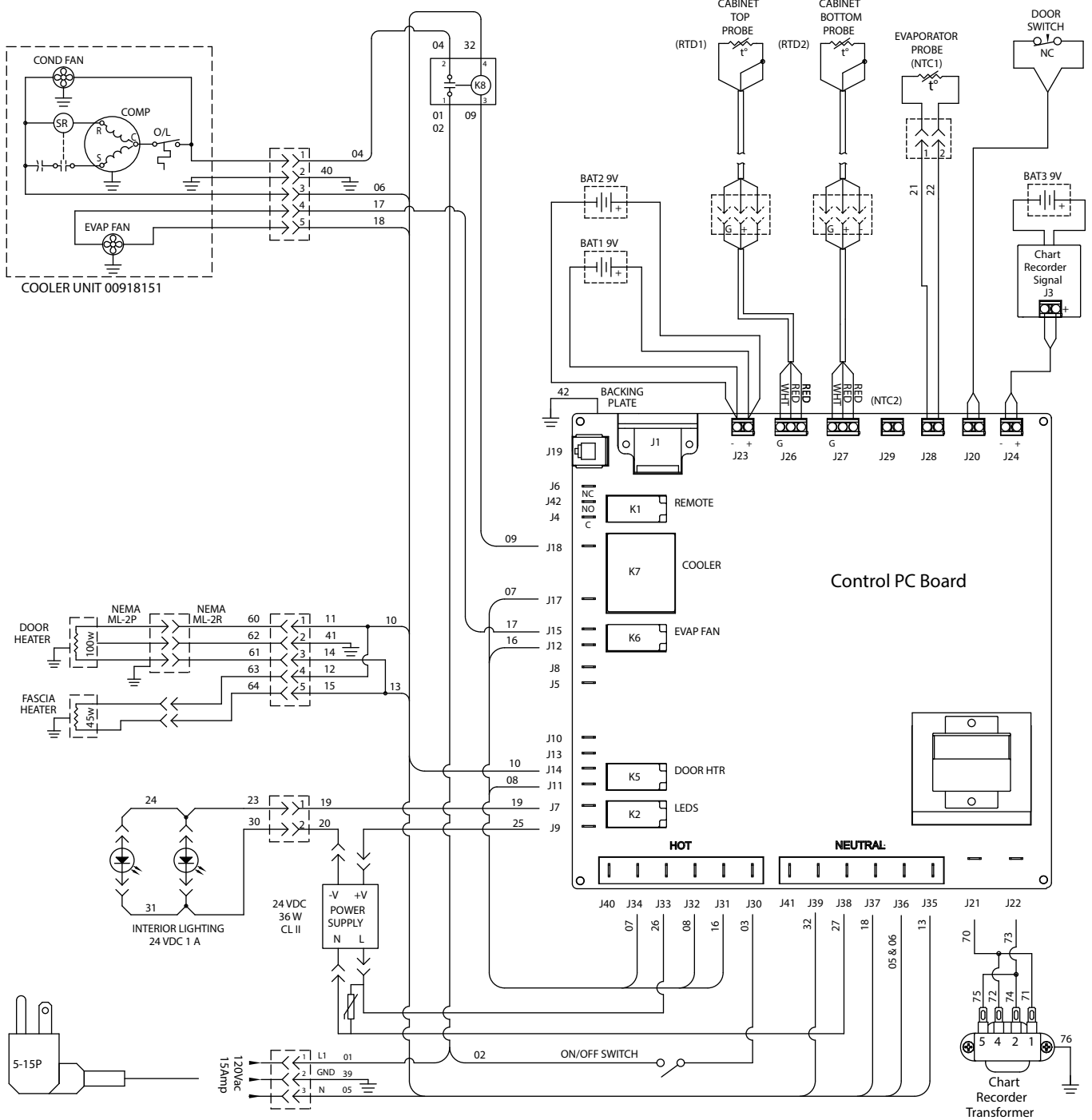

Component Details

Compressor amperage

RLA - 7.6A, LRA - 44.5A

Compressor winding resistances

C-S (5.2Ω - 5.9Ω), C-R (1.2Ω - 1.3Ω)


Evaporator fan motor amperage

FLA - 0.22A


Condenser fan motor amperage

FLA - 0.46A

Wiring diagram (Serial Numbers D61742— E15526)

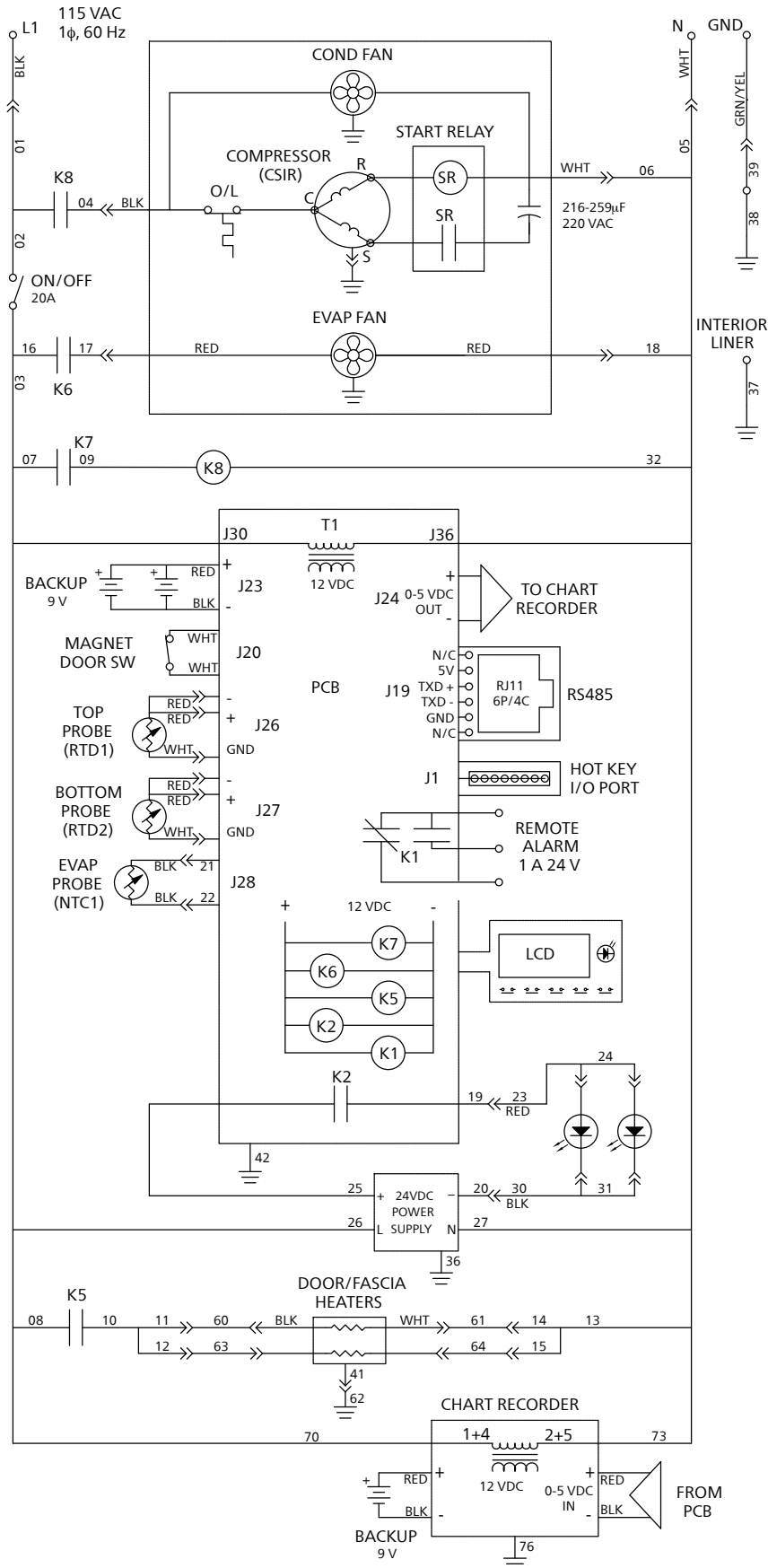

Wiring diagram (Serial Numbers D61742— E15526)

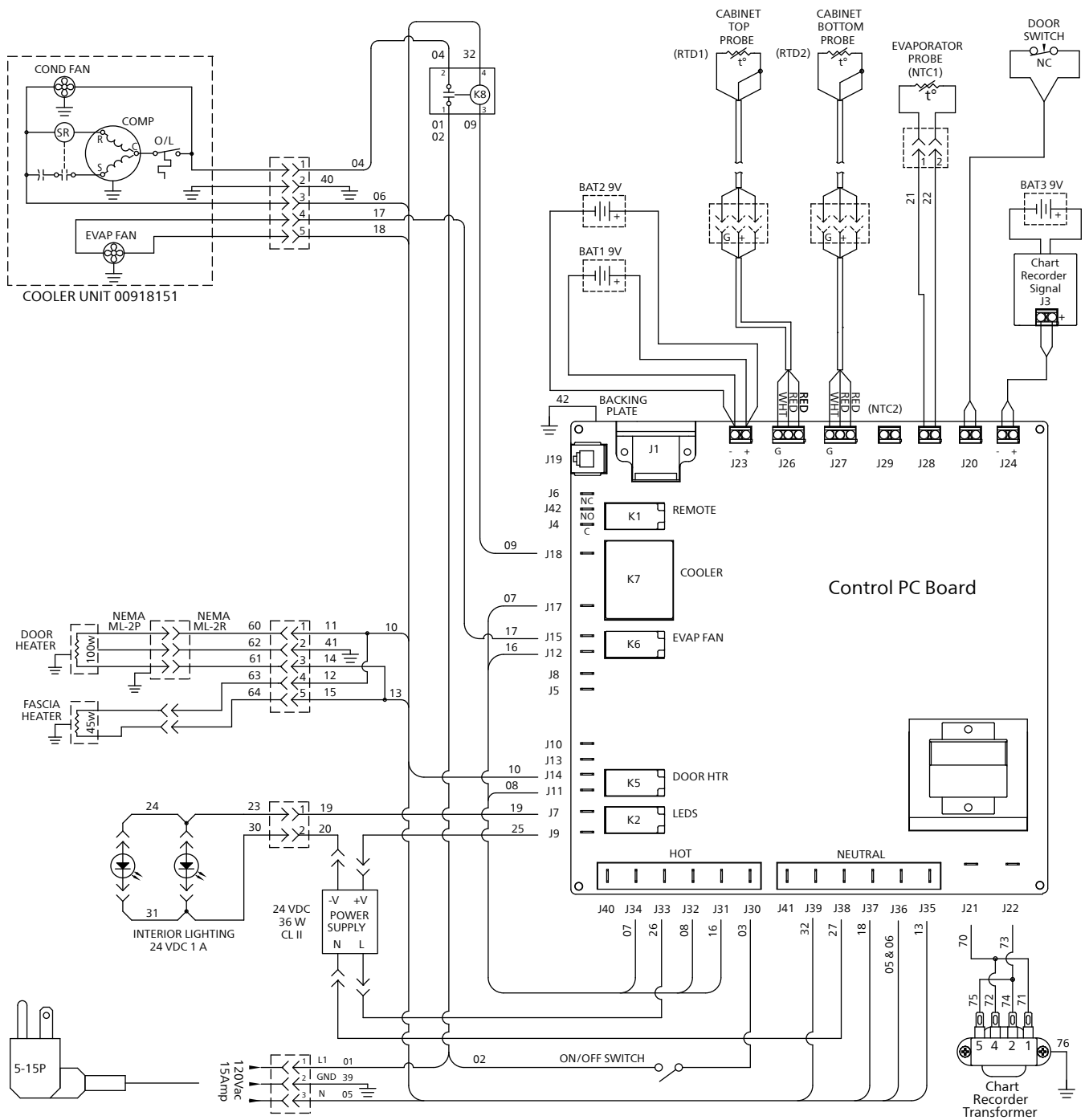

Component Details

Compressor amperage

RLA - 7.6A, LRA - 44.5A

Compressor winding resistances

C-S (5.2Ω - 5.9Ω), C-R (1.2Ω - 1.3Ω)


Evaporator fan motor amperage

FLA - 0.22A


Condenser fan motor amperage

FLA - 0.46A

Wiring diagram (Before Serial Number D61742)

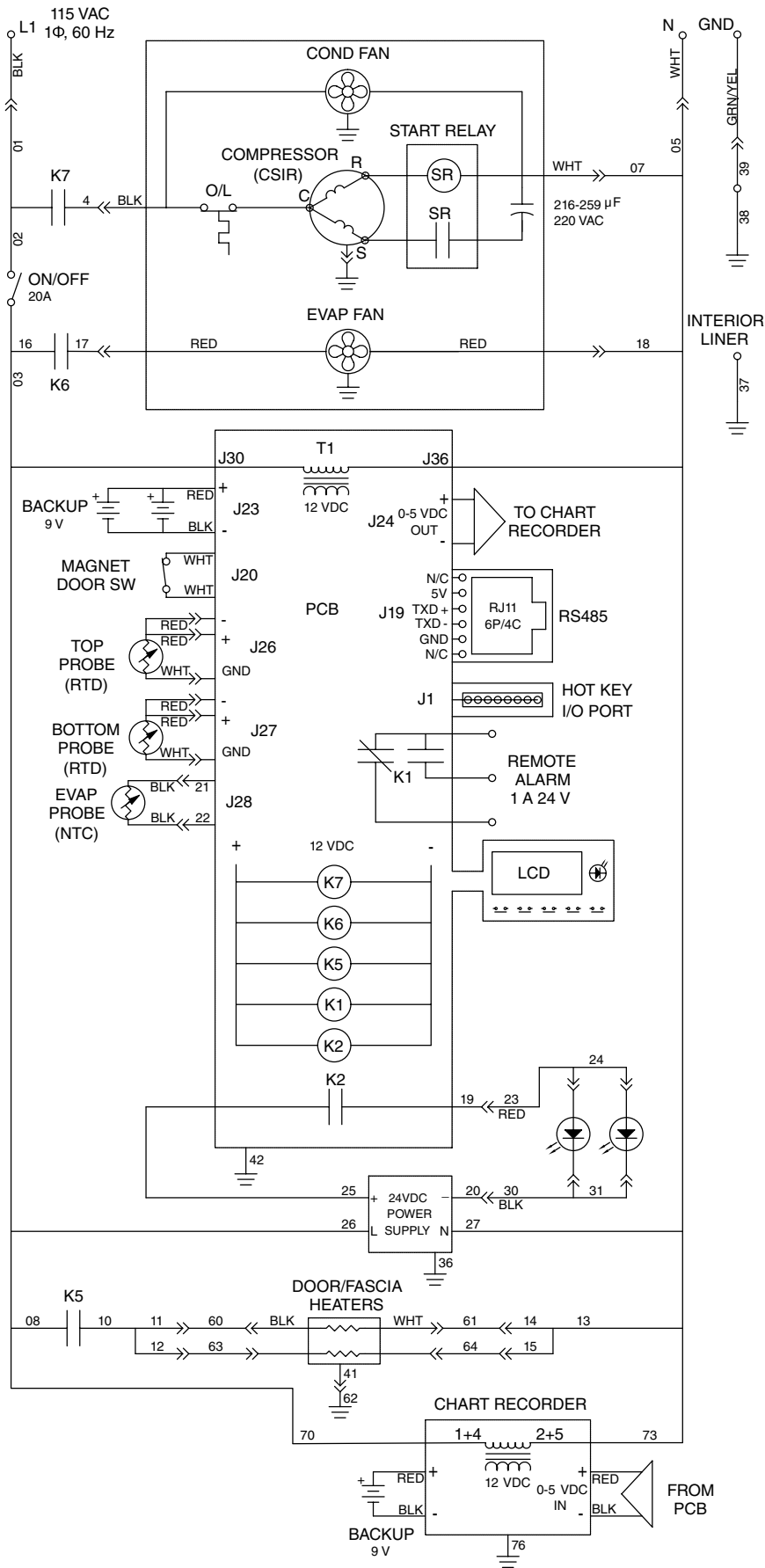

Wiring diagram (Before Serial Number D61742)

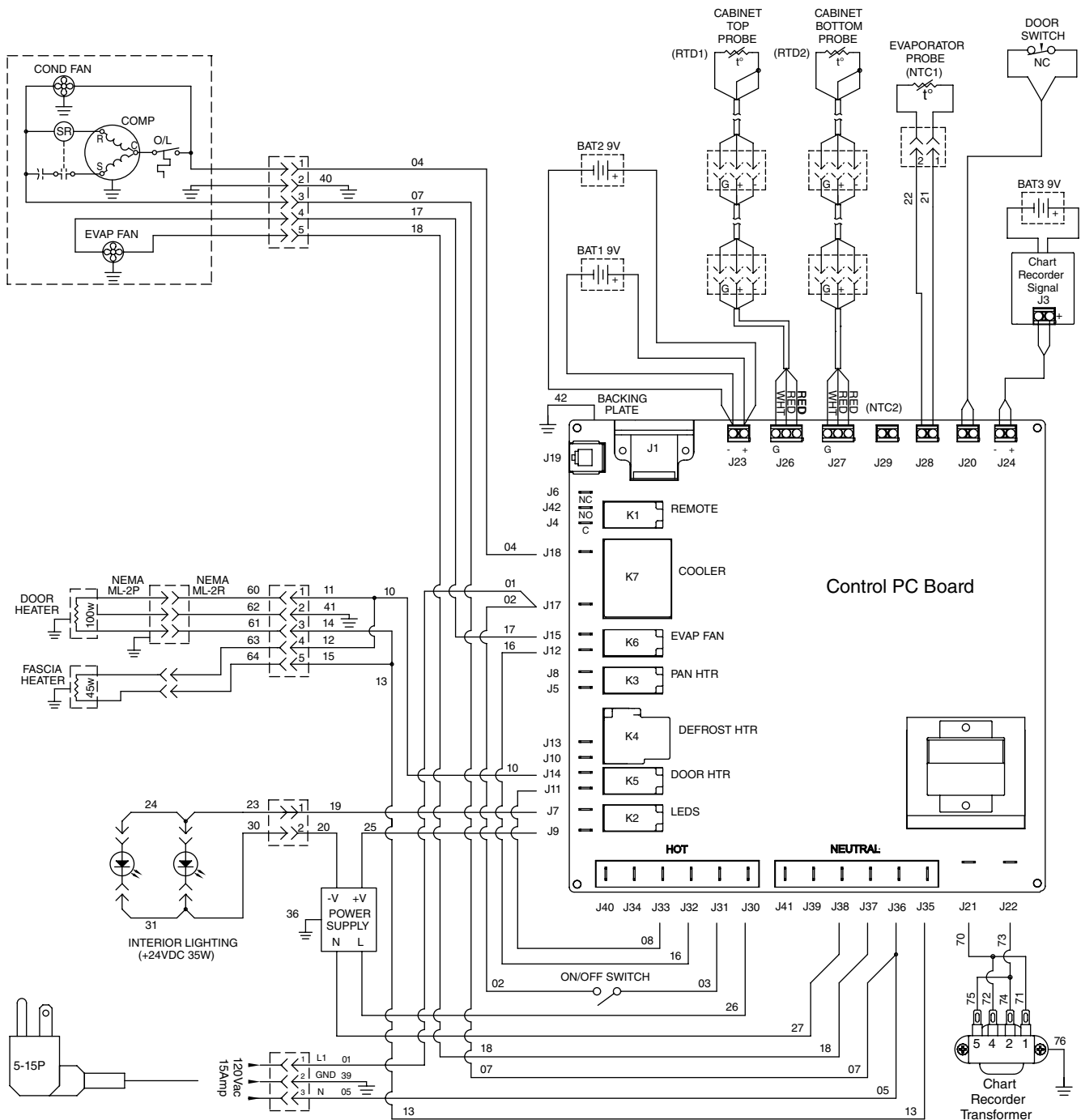

Component Details

- Compressor amperage
RLA - 7.6A, LRA - 44.5A
- Compressor winding resistances
C-S (5.2Ω - 5.9Ω), C-R (1.2Ω - 1.3Ω)
- Evaporator fan motor amperage
FLA - 0.22A
- Condenser fan motor amperage
FLA - 0.46A

Troubleshooting

Before calling for service


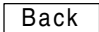







1. Check that unit is plugged in.
2. Test outlet with another appliance to verify power.

Problem	Indicators	Corrective Action
Refrigerator does not operate (no components run).	1. Power switch faulty or in OFF position; loose connection.	1. Turn power switch to ON position; check switch and connections.
	2. Refrigerator not plugged in.	2. Connect plug.
	3. No power at outlet.	3. Restore power.
	4. Temp controller not energizing components.	4. Check controller contact terminals for power. Replace controller if needed.
	5. Probe not sensing cut-in temperature.	5. Replace controller and/or probe.
Compressor does not run.	1. Thermal overload open or defective.	1. Allow to cool or replace.
	2. Capacitor and/or relay defective.	2. Replace as required.
	3. Compressor defective.	3. Replace compressor.
Evaporator fan motor does not run.	1. Defective fan motor.	1. Replace fan motor
Refrigerator does not shutoff.	1. Controller not sensing cut of temperature.	1. Replace controller and/or evaporator (NTC1) probe.
	2. Controller keeping refrigeration system energized.	2. Replace controller.
Refrigerator does not maintain temperature (all components run).	1. Condenser or evaporator coil needs cleaning.	1. Clean coils as needed.
	2. Door not closing.	2. Adjust door closure per page 5.
	3. Excessively high ambient.	3. Maximum recommended ambient is 100 F.
	4. Refrigerant leak.	4. Locate and repair leak.
	5. Incorrect refrigerant charge.	5. Recover, evacuate and weigh in correct charge.
	6. Plugged capillary tube.	6. Replace capillary tube.
	7. Inefficient compressor.	7. Consult technical service.
	8. Faulty door gasket.	8. Replace door gasket

If problems persist after following this basic troubleshooting guide, call Follett's technical service group at (877) 612-5086 or +1 (610) 252-7301.

Appendix

Control key functions


 MENU	Enter main menu	 Back	Return to prior		Buttons to select screen options
 ▲	Scroll up	 ENTER	Select		Cabinet light ON/OFF
 ▼	Scroll down	 MUTE	Mutes alarm		Alarm light

Menu Navigation table

Main Menu	Sub menu	Default setting	Range	Notes
If password is ON, enter 4 digit pin		3843	0000-9999	Adjust with up/down arrow, next moves to next digit, press enter to set
Alarm settings	High alarm set	40 C (104 F)	40 to -40 C (104 to -40F)	Default is set high to not alarm on initial cool down
	Low alarm set	1.0 C (33.8 F)	40 to -40 C (104 to -40F)	
Display settings	Show bottom temp	NO	YES/NO	
	Display inverse	YES	YES/NO	Reverse light lettering with dark background
	Degrees F/C	C	F or C	Displays temperatures in F or C
Min/Max temp				Records min/max temperature and time of top probe. Up/down arrow displays min/max. Enter to reset.
Batt = _V OK				Shows status of (2) 9V batteries.
Refrigeration controls (press and hold down-arrow and ENTER for 3 seconds)	Calibration	Cal top probe		Adjust with up/down arrow, press enter to set
		Cal bottom probe		Adjust with up/down arrow, press enter to set
	Refrig. setpoint	0.5 C (32.9 F)	40 to -40 C (104 to -40F)	
	Differential	5.0 C (9.0 F)	3.0 to 6 C (5.4 to 10.8 F)	
	High setpoint limit	5.0 C (41.0 F)	40 to -40 C (104 to -40F)	
	Low setpoint limit	-1.5 C (29.3 F)	40 to -40 C (104 to -40F)	
	Door heater cycle	0%	0 to 100%	Heater on percent (50% for heated door option)
	On time	4 min	2-14 min	Compressor on time
	Off time	8 min	2-14 min	Compressor off time
	Temp sensors			Displays probe temperatures
Password ON/OFF		OFF	ON/OFF	
PIN		3843	0000-9999	Any four digit PIN can be stored beside the permanent factory default. Adjust with up/down arrow, next moves to next digit, press enter to set.
Set Time/Date				Adjust with up/down arrow, next moves to next digit, press enter to set
Set alarm volume		High	High/Med/Low	Buzzer volume
Alarm silence		ON	ON/OFF	Audible alarm


Replacement parts

Unit cooler

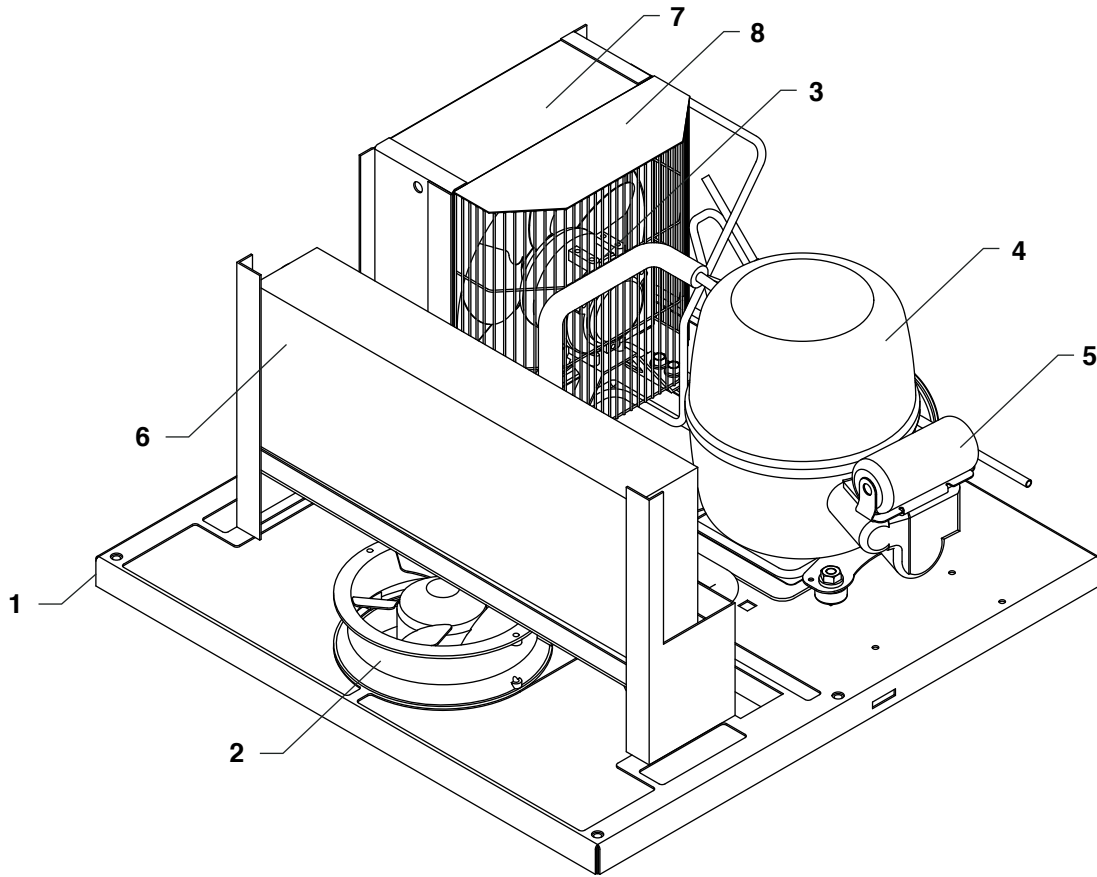


Reference	Description	Part #
Not shown	Gasket, unit cooler	00912980
Not shown	Probe, evap, NTC	00940148
1	Unit cooler with mounting bolts and gasket	00939785
2	Evaporator fan motor with blade	00945311
3	Condenser fan motor with blade	01179704
Not shown	Cap tube and drier	00945378
4	Compressor with start capacitor, relay and overload	01096015
5	Start capacitor	01077734
Not shown	Start relay	00945360
Not shown	Compressor overload	00945386
6	Evaporator	00945766
7	Condenser	00945774
8	Fan guard, condenser	00945782
Not shown	Cord and plug kit	00939900
Not shown	Terminal cover, compressor	01223726

Cabinet

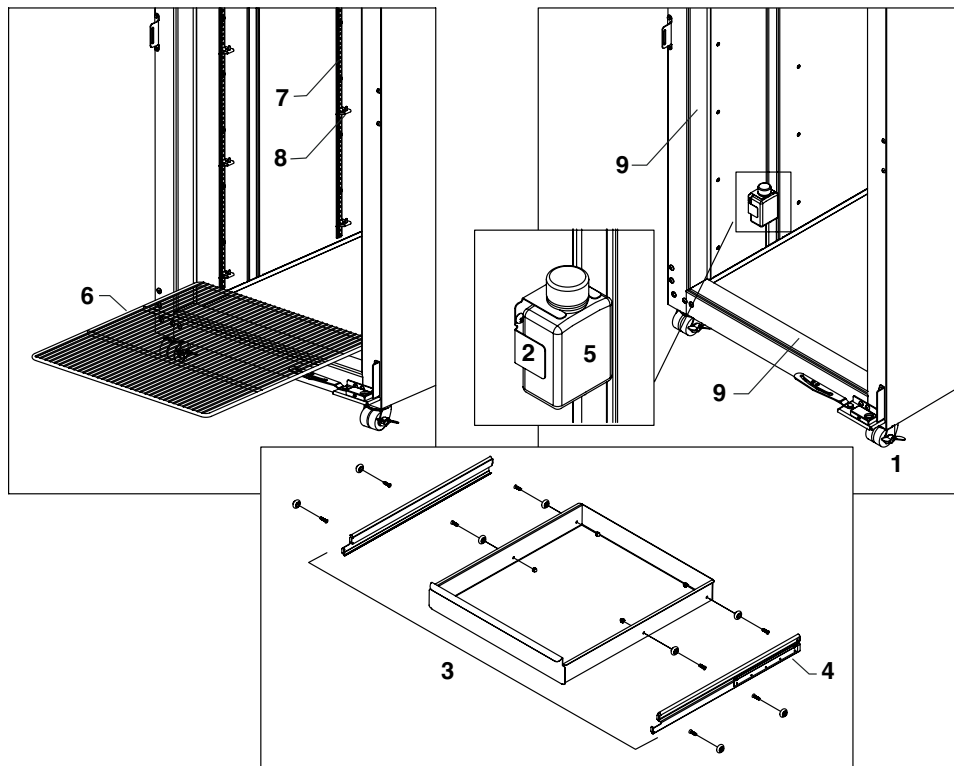


Reference	Description	Part #
1	Caster, swivel with brake	00187674
2	Bracket, bottle, 125ML	00171132
3	Drawer, assembly, REF25	00939652
3	Drawer, assembly, REF20	00939660
4	Slide, drawer, REF25 (set of 2)	00193045
4	Slide, drawer, REF20 (set of 2)	00188367
5	Bottle with gasket	01015817
Not shown	Basket, assembly, REF25	00939710
Not shown	Basket, assembly, REF20	00939728
6	Shelf, REF25 (with supports)	01015841
6	Shelf, REF20 (with supports)	01015858
7	Pilaster	00901173
Not shown	Pilaster, screws	00902213
8	Support, shelf	00156240
Not shown	Air panel	00901561
Not shown	Probe, RTD, top	00918839
Not shown	Probe, RTD, NIST calibrated, top	01101096
Not shown	Probe, RTD, bottom	00988055
Not shown	Probe, RTD, NIST calibrated, bottom	01101104
Not shown	Plug, RTD probe	00919274
Not shown	Panel jack, female	00918847
9	Cover, thermal break (4 pieces)	00922997
Not shown	Channel, probe	01004159
Not shown	Bushing, probe channel (LB and PH require 2, BB requires 3)	00111104
Not shown	LED, light strip	00980912
Not shown	Wire clip, probe	00919910
Not shown	Glycerine (1 pint)	00959296
Not shown	LEDs (2), MOV and instructions	01049709

Facade

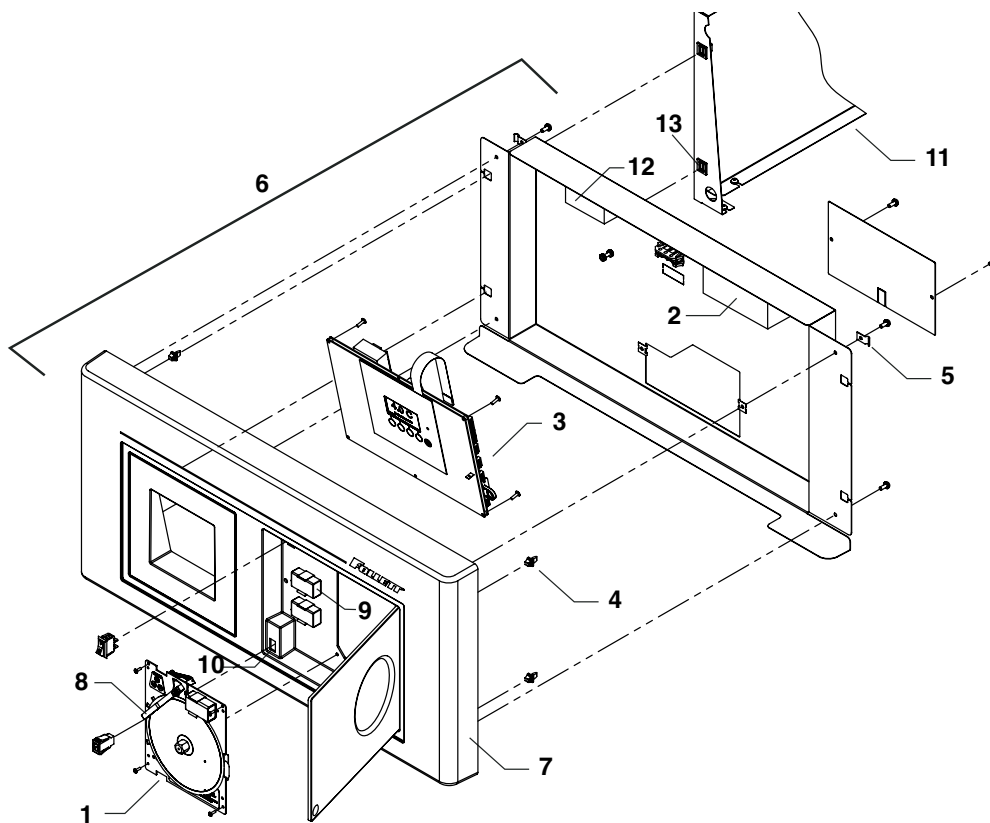

Reference	Description	Part #
1	Chart recorder	00906198
2	LED power supply	00939801
3	Controller (for serial numbers above D61742)	01144138 or 01144146
3	Controller (for serial numbers below D61742)	01144138 or 01144146
Not shown	Kit, door switch	00939959
4	Male, bayonet (strike, snap)	00179549
5	Facade, locking clip	00931501
6	Complete facade assembly without chart recorder	01144138
6	Complete facade assembly with chart recorder	01144146
7	Facade, cover, with chart recorder	00940130
Not shown	Facade, cover, without chart recorder	00940155
8	Pens, chart recorder, red (box of 6)	00918813
Not shown	Charts (box of 50) in degrees C	00918821
Not shown	Charts (box of 50) in degrees F	00967729
9	Battery, 9V	00112177
10	Latch, touch	00901165
11	Shell, upper side REF25	00940098
11	Shell, upper side REF20	00925933
Not shown	Top ventilated cover	00991760
12	Control relay, compressor (for serial numbers above D61742)	00980482
Not shown	Bottom probe, resistor	01041342
Not shown	MOV assembly	01042258
Not shown	Back support	00991695
13	Female, fastener	00179556

Door assembly

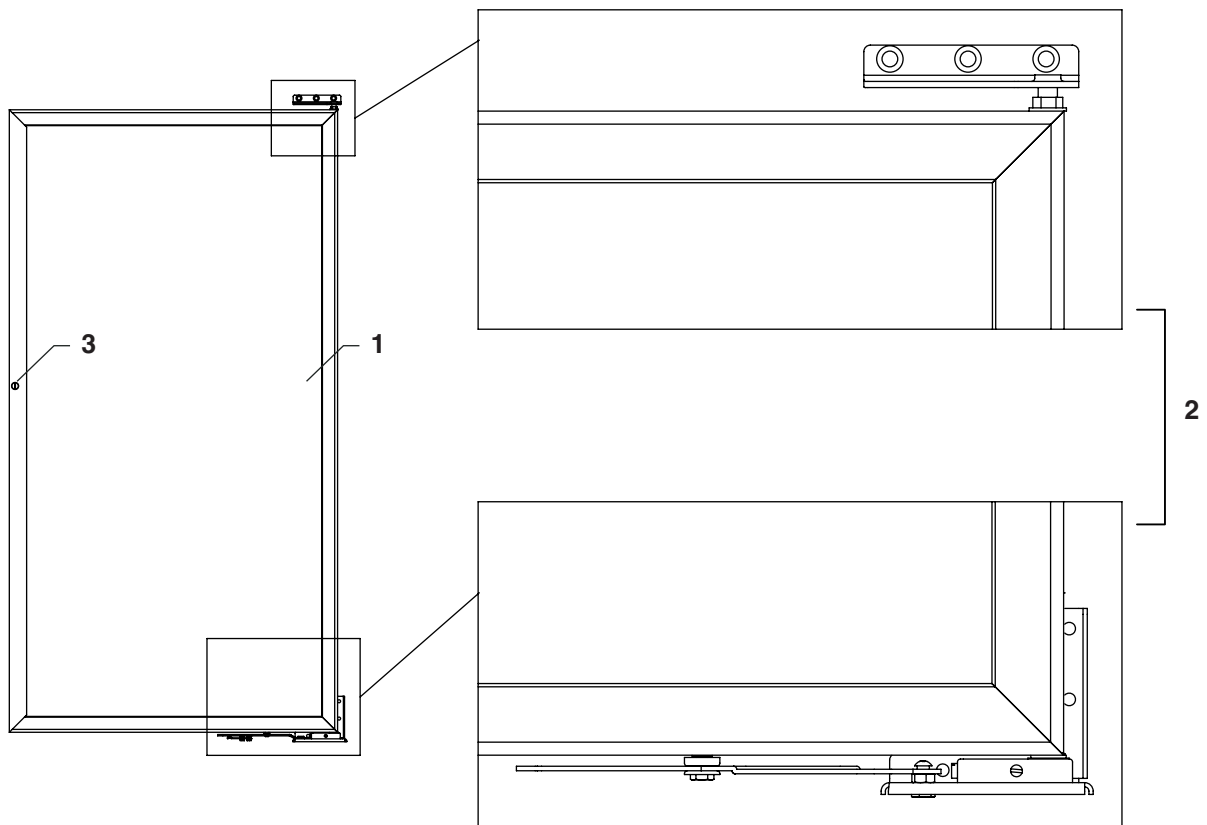

Reference	Description	Part #
Not shown	Door gasket	00945345
Not shown	Left-hand, non-heated door	00939876
1	Right-hand, non-heated door	00939884
2	Right-hand hinges and hardware with black colored hold open bracket	00996579
Not shown	Left-hand hinges and hardware with black colored hold open bracket	00996587
Not shown	Left-hand hinges and hardware with steel colored hold open bracket	00939918
Not shown	Right-hand hinges and hardware with steel colored hold open bracket	00939926
Not shown	Door magnet kit	00939967
Not shown	Bracket, strike	00939983
Not shown	Left-hand heated door	00939991
Not shown	Right-hand heated door	00940007
Not shown	Key, door	00945436
3	Lock, door	00945444
Not shown	Soft port kit	00940122
Not shown	Wall bracket kit, seismic anchoring	00927194
Not shown	Kit, Pyxis bracket	00927202
Not shown	Kit, Omnicell	00966432

Warranty Registration and Equipment Evaluation

Thank you for purchasing Follett® equipment. We hope you find that our equipment meets or exceeds your expectations, as our goal is to deliver high value products and services that earn your complete satisfaction!

Please review the enclosed installation and operations manual. It is important that the installation be performed to factory specifications, so your equipment operates to its maximum efficiency.

Follett LLC will not be liable for any consequential damages, expenses, connecting or disconnecting charges or any losses resulting from a defect of the machine.

For full warranty details, visit our website www.follettice.com/productwarranties.

Warranty registration and equipment evaluation is important to help us keep track of our equipment and to record the machine's performance. We request that you register Follett equipment warranties on our website www.follettice.com/support and choose Warranty Registration and Equipment Evaluation. It's simple to do; please take a moment to register today. There is also space on the form to provide us with comments and feedback. Please let us know about your experience so we can capture it for our continuous improvement efforts.

We pride ourselves on producing outstanding equipment and we work hard to back it up with outstanding customer and technical support. Please let us know what else we can do to assist you. We would be happy to answer your questions.

Registro de Garantía y Evaluación del Equipo

Gracias por haber elegido este producto Follett®. Esperamos que nuestro equipo cumpla o supere sus expectativas porque es nuestro objetivo ofrecer productos y servicios de gran valor que se ganen su plena confianza.

Le rogamos consulte el manual de instalación y de instrucciones adjunto, ya que es muy importante que la instalación se realice según las especificaciones de fábrica para que el equipo funcione a su máxima eficiencia.

Follett LLC no se hace responsable de los daños indirectos, costos, gastos por conexión y desconexión o pérdidas por causa de defecto de la máquina.

Si desea una información más completa sobre la garantía, visite nuestro sitio web www.follettice.com/productwarranties.

Las operaciones de registro de la garantía y evaluación del equipo son importantes para que podamos realizar un seguimiento de nuestro equipo y registrar el rendimiento de la maquinaria. Por favor, registre las garantías del equipo Follett en nuestro sitio web www.follettice.com/support y seleccione Registro de la Garantía y Evaluación del Equipo. Es muy sencillo, solo le llevará un momento realizar hoy mismo el registro. En el formulario incluimos un espacio en blanco para sus comentarios y opiniones. Infórmenos sobre su experiencia para que podamos incorporarla a nuestros continuos esfuerzos de mejora.

Nos enorgullecemos de producir un equipo excepcional y trabajamos duro para respaldarlo con un soporte técnico y un servicio de atención al cliente de primera. Le rogamos nos indique qué más podemos hacer para ayudarlo. Estaremos encantados de responder a sus dudas.

Enregistrement de la garantie et évaluation de l'équipement

Merci d'avoir acheté un équipement Follett®. Notre objectif étant d'offrir des produits et des services de grande valeur vous satisfaisant pleinement, nous espérons que celui-ci satisfera, voire dépassera, vos attentes !

Veuillez consulter le manuel d'installation et d'exploitation. Il est important que l'installation soit réalisée conformément aux spécifications de l'usine, de sorte que votre équipement fonctionne à son rendement maximum.

Follett LLC n'est pas responsable de tout dommage consécutif, de toute dépense, de tout frais de raccordement ou de déconnexion, ni de toute perte liée à un défaut de la machine.

Pour lire la garantie dans son ensemble, visitez notre site Internet www.follettice.com/productwarranties.

L'enregistrement de la garantie et l'évaluation de l'équipement sont importants pour nous aider à suivre notre équipement et pour enregistrer les performances de la machine. Nous vous demandons donc d'enregistrer la garantie de votre équipement Follett sur notre site Internet, www.follettice.com/support, dans la section Warranty Registration and Equipment Evaluation. Cette opération est simple ; veuillez prendre un moment pour la réaliser aujourd'hui.

Le formulaire contient également un espace pour nous faire parvenir vos commentaires et un retour d'informations. Veuillez nous faire part de votre expérience pour que nous puissions prendre appui dessus pour poursuivre nos efforts constants d'amélioration.

Nous sommes fiers de produire des équipements exceptionnels et nous travaillons avec acharnement pour y associer une assistance à la clientèle et technique exceptionnelle. N'hésitez pas à nous indiquer dans quelle mesure nous pouvons vous aider. Nous serions ravis de répondre à vos questions.